PC controlled stirring

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The screenshot displays the software interface for process control, divided into three main sections:

- Grafik (Graph):** Located on the left, it features a coordinate system with a vertical axis from 0 to 120 and a horizontal axis labeled 'Time (minutes)' from 00:00 to 00:00. A 'Menüleiste' (menu bar) is positioned above the graph.
- Prozessparameter (Process parameters):** Located on the right, it shows a 'Process loading' section with a device image and a 'Parameter settings' panel. The settings include:
 - Time: 00:00
 - Start: 00:00:00
 - Start speed: 2000 rpm
 - Actual speed: 10 rpm
 - Torque: 100%
- Benutzerprofile (User profiles):** Located at the bottom, it contains a 'Profile management' table with columns for 'Name', 'Speed', and 'Load Weight'. Below the table is a large empty grid area.

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Translation of the original instructions
Page 32 – 57

Content

Introduction

About this document	32
Typographic conventions	32
Copyright protection	32

General notes

Basic product information	33
Residual risk	33
Intended use	33
Reasonably foreseeable misuse	33

Safety

General safety information	34
Operational safety	34
Other regulations	34

Installation

Installation instructions	35
General notes	35
System requirements	35
Install the USB driver	36
Install Hei-CONTROL	38
Update/adjust/restore Hei-CONTROL	42

Commissioning

Commissioning	45
Start the software	45
Start device(s)	45
Communication	45

Operation

User interface	46
Home screen	46
Process parameters screen	47
Process view	48
Control elements	49
Graphics area	50
Process parameters area	52
Monitor devices	53
Control devices	53
User profiles area	54

Troubleshooting

Troubleshooting	56
Contact information Heidolph international	57
Warranty Statement	57

About this document

These operating instructions describe all functions and operation of the Hei-CONTROL software for controlling Hei-TORQUE overhead stirrers as well as Hei-CONNECT and Hei-END magnetic stirrers manufactured and distributed by Heidolph instruments GmbH & Co. KG. The operating instructions manual is an integral part of the delivery!

Typographic conventions

Standardized symbols, highlighting elements, and signal words are used in this document to identify warnings, cautions, important information, and special text contents.

Symbol	Signal word / explanatory note
	<p>Warning symbols in combination with a signal word indicate dangers:</p> <p>DANGER Indicates a hazardous situation which, if not avoided, will result in death or serious injury.</p> <p>WARNING Indicates a possible hazardous situation which, if not avoided, may result in death or serious injury.</p> <p>CAUTION Indicates a hazardous situation which, if not avoided, could result in minor or moderate injury or material/environmental damage.</p>
	<p>Prohibition signs indicate actions or situations that are to be omitted or avoided.</p> <p>Failure to comply may result in personal injury and / or damage to property.</p>
	<p>Mandatory signs are used to indicate important information regarding the product handling.</p> <p>This information is used to ensure operational safety and to maintain the value of the product.</p>
[GUI]	<p>Parameter designations, display texts, and device labels are highlighted in text and tables in a typographic manner to facilitate the assignment on the device.</p>
→	<p>The arrow symbol indicates instructions to be followed in order to ensure the operational safety when handling the product.</p>

Copyright protection

This publication is protected by copyright and intended for internal use by the purchaser of the product only.

No part of this publication may be transmitted or reproduced in any form, by any means, without the prior written consent of the copyright owner Heidolph Instruments GmbH & Co. KG. Any violation is subject to compensation for damages.

Basic product information

Residual risk

The product was designed in accordance with the latest technical standards at the time of development and the recognized safety regulations.

The product described does not pose any direct risks.

When using the product in accordance with the regulations for proper use, all operating instructions and safety instructions for the controlled overhead or magnetic stirrers must be observed in order to avoid hazards and to ensure general operating and occupational safety (see the manufacturer's operating instructions)!

Intended use

The Hei-CONTROL software was designed exclusively for controlling Hei-TORQUE overhead stirrers as well as Hei-CONNECT and Hei-END magnetic stirrers manufactured and distributed by Heidolph Instruments GmbH & Co. KG.

Any other use of this software is not considered as intended!

Reasonably foreseeable misuse

Compliance with and implementation of all relevant guidelines and safety measures for the respective field of application is within the sole responsibility of the operator.

All risks resulting from improper use of the product are solely borne by the operator.

The product may exclusively be operated by authorized and instructed personnel. Training and qualification of the operating personnel as well as ensuring that the product is operated with responsibility are the sole responsibility of the operator!

General safety information

- Before installing and using the software, familiarize yourself with all the safety regulations and guidelines for occupational safety applicable at the place of use and observe them at all times.
- Before using the software, familiarize yourself with all regulations and guidelines applicable for the controlled device(s) and observe them at all times.
- If there is missing or misleading information on the product or on occupational safety, contact the responsible safety specialist or our technical service.
- Only use the software in accordance with the regulations for intended use (see section „Intended use“ on page 33).

Operational safety

- Do not make any arbitrary changes to the software, do not reprogram the software by yourself!
- In the event of an error, immediately disconnect all controlled devices (close the program on the PC, disconnect the cable connection between PC and device), switch off the controlled devices and secure them against unintentional restart.
- Before resuming operation, ensure that any software malfunctions are resolved.
- Contact our technical service if troubleshooting is not immediately possible (see section „Contact information Heidolph international“ on page 57).

Noncompliance will invalidate any warranty against Heidolph Instruments.

The operator is exclusively liable for all damages resulting from unauthorized changes to the software or from disregard of the safety and danger instructions or the instructions for use of the manufacturer!

Other regulations

In addition to the notes and instructions in this document, observe all other applicable regulations such as laboratory and workplace guidelines, hazardous substances ordinances, recognized rules of safety engineering and occupational medicine as well as particular local regulations!

Installation instructions

General notes

Install the necessary USB driver and software before connecting the PC to the device(s)!

In case of an update, the existing program version must first be uninstalled using the installation routine!

It is not possible to update the control software by overwriting it.

System requirements

The Hei-CONTROL software can be installed on any PC that meets the following system requirements:

System requirement	Minimum	Recommended
Operating system	Windows®7 or later, 32-bit	Windows®7 or later, 64-bit
Memory	2 GB	4 GB
Processor	2.5 GHz dual-core processor	2.5 GHz 4-core processor
Interfaces	1×USB or 1×RS232	depending on the number of units controlled
Display	6,500 colors	16 million colors
Screen resolution	1.400 × 1.050 pixels	1.680 × 1.050 pixels

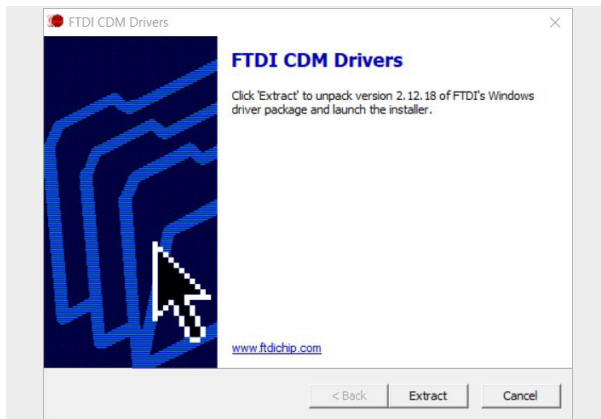
Install the USB driver

Hei-TORQUE laboratory stirrers are equipped with a USB type interface which requires a special USB driver to be installed on the controlling PC.


In case of missing system authorizations (admin rights), please contact the responsible system administrator!

- Connect the supplied USB stick to an USB port of your PC.
- Open the Windows® Explorer and access the **Hei-Control_Vx** folder stored on the stick (x = version-dependent).
- Open the subfolder **usb driver**.
- Start the driver installation by double-clicking the **CDM vx WHQL Certified.exe** file (x = version-dependent):


- Click **Extract** to unpack the driver package.

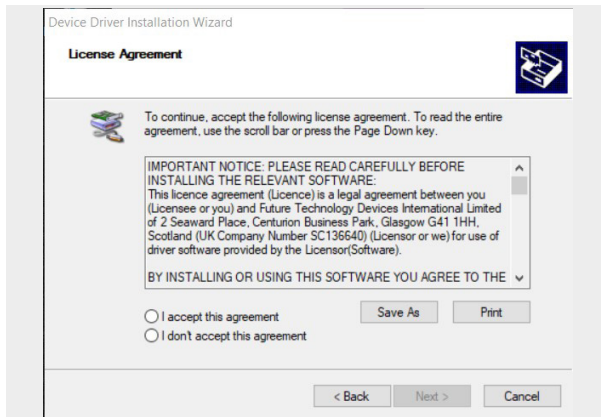- After extraction, the installation wizard opens. Click **Next** to open the **License Agreement** dialog box:

- Read and confirm the license agreement to proceed with the driver installation: Select **I accept this agreement** and continue the driver installation with **Next**.
- You have the option of storing the license agreement in a location of your choice (**Save As**) or print from a connected printer (**Print**).
 - The installation routine can be canceled at any time with **Cancel**.
- When the driver installation is complete, the following dialog box appears:

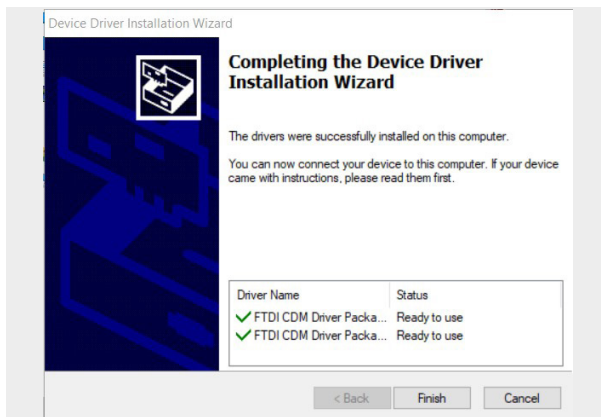- Close the installation routine by clicking **Finish**.

Install Hei-CONTROL

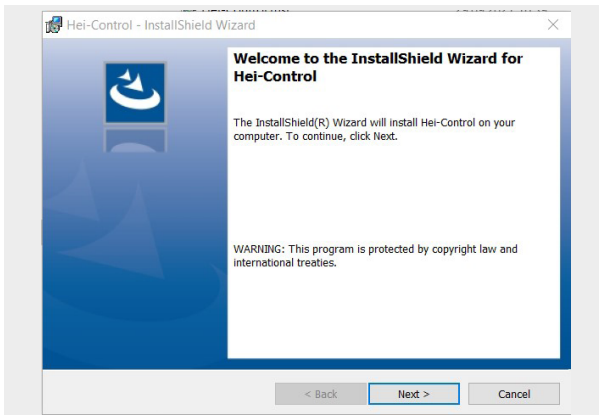
This section describes the procedure for installing the Hei-CONTROL control software for the first time.

The software installation can only be performed with the appropriate authorization on your PC (Windows® option **Run as administrator!**)


In case of missing system authorizations (admin rights), please contact the responsible system administrator!

- Connect the supplied USB stick to an USB port of your PC.
- Open the Windows® Explorer and access the **Hei-Control_Vx** folder stored on the stick (x = version-dependent).
- Start the installation by double-clicking on the **setup.exe** file:

- On the start screen, confirm with **Next** to continue.

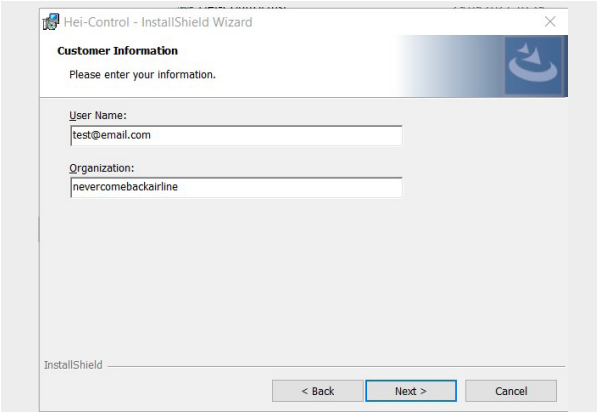When the software is reinstalled, the **Customer Information** screen appears:

Hei-Control - InstallShield Wizard

Customer Information

Please enter your information.

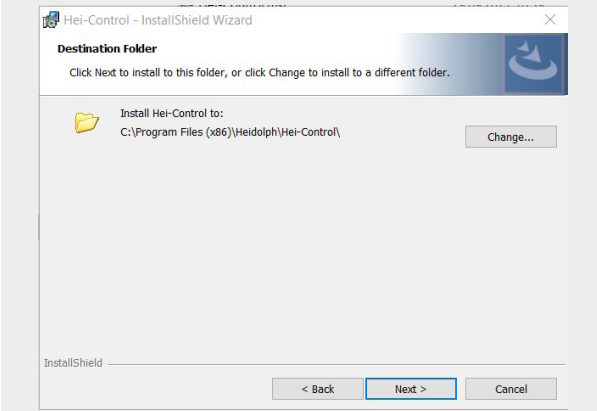
User Name:
test@email.com

Organization:
nevercomebackairline

InstallShield

< Back Next > Cancel

- Enter your user information in the **User Name:** and **Organization:** fields.
- Confirm with **Next** to open the following **Destination Folder** screen.
- On the **Destination Folder** screen, specify a destination folder.

Hei-Control - InstallShield Wizard

Destination Folder

Click Next to install to this folder, or click Change to install to a different folder.

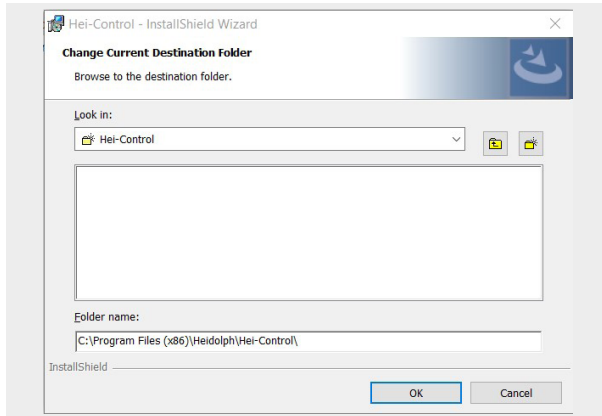
Install Hei-Control to:
C:\Program Files (x86)\Heidolph\Hei-Control\

Change...


InstallShield

< Back Next > Cancel


- Click **Next** to accept the default destination folder.
- or**
- Click **Change** to specify a user-specific destination folder:

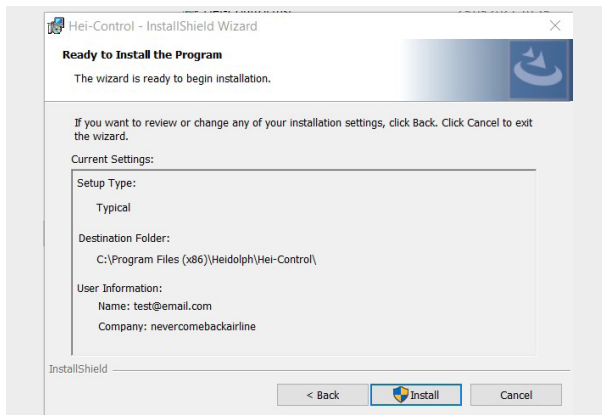- On the **Change Current Destination Folder** screen, specify the destination folder you want to use.
- Confirm with **OK** to return to the destination **Destination Folder** screen.
- On the **Destination Folder** screen, confirm the new destination folder by clicking **Next**.
- Confirm the summary by clicking **Install** to start the installation process:

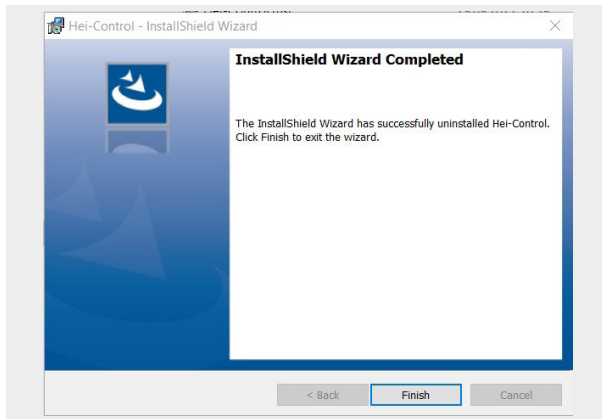- Once all software components have been successfully installed, the installation routine completion window appears:

- Click **Finish** to exit the wizard. The software is ready for use without restarting the PC.

Start the software by double-clicking on the program icon on the desktop of your PC or via the two icons (**Hei-Control.exe**, **MR-Control.exe**) in the program folder.

During installation, these operating instructions are also stored as a PDF file in the program folder.

Update/modify/restore Hei-CONTROL

This section describes how to update, customize, or repair the Hei-CONTROL software.

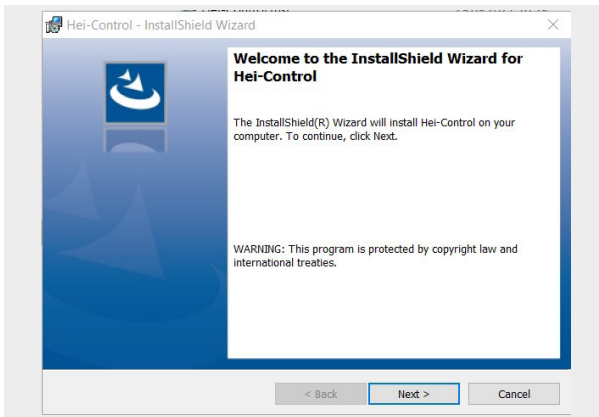
The existing software version must be uninstalled before updating to a newer version. Use the **Remove** option of the installation wizard or the standard uninstall routine of Windows®!

Existing profiles are retained in the program folder when the software is uninstalled.


The software can only be updated, customized, or removed with the appropriate authorization on your PC (Windows® option **Run as administrator**)!

In case of missing system authorizations (admin rights), please contact the responsible system administrator!

- Start the installation routine by double-clicking on the **setup.exe** file stored in the program folder on the hard disk of the PC:

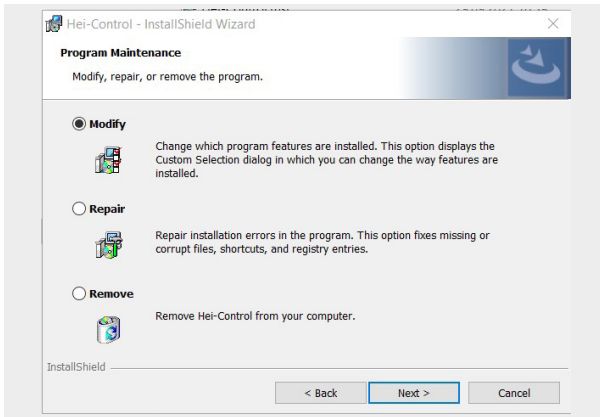- On the start screen, confirm with **Next** to continue. The **Program Maintenance** screen opens:

→ The following options are available:

Modify

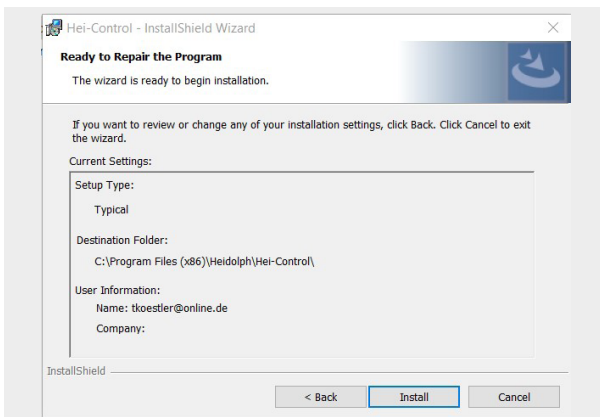
The **Modify** option is not supported in the current version.

→ Return to the **Program Maintenance** screen by clicking **Back**.

Repair

→ To repair the program, select **Repair**.

→ Click **Next** to open the **Ready to Repair the Program** screen:


→ Confirm the summary with **Install** to reinstall the components of the current version:


→ Return to the **Program Maintenance** screen by clicking **Back**.

Remove

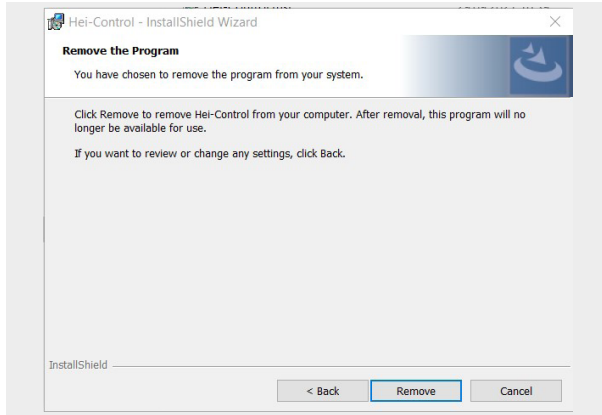- To uninstall the current program version before updating, select **Remove**.
- Press **Next** to continue.
- In the **Remove the Program** screen, confirm by clicking **Remove** to remove the existing program version from your PC:

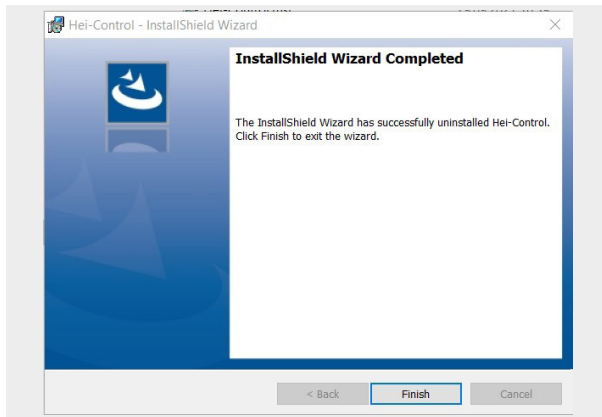- Once all software components have been successfully removed, the installation routine completion window appears:

- Confirm with **Finish** and follow the instructions in section „Install Hei-CONTROL“ on page 38 to install the new version of the program.

Commissioning

After installing the driver and software, connect the device(s) to the PC you are using. The available interfaces and supplied cables vary depending on the device and version. Observe the operating instructions supplied with the individual device (PC, laboratory stirrer, magnetic stirrer)!

Any necessary adapters (USB to RS232) are not included in the scope of delivery and must be obtained from a specialist dealer.

Start the software

CAUTION

Possible personal injury/property damage and/or production loss due to unexpected start-up of the device

- Before starting the software, make sure that all connected devices are turned off.
 - Before starting the software, check all cables and connectors for damages and make sure that they are securely attached.
 - Observe the operating instructions(s) supplied with the individual device and follow all safety instructions and work instructions contained therein!
-

Start the control software by double-clicking on the program icon on the desktop of your PC or via the two icons (**Hei-Control.exe**, **MR-Control.exe**) in the program folder.

Make all necessary settings in the user interface, see section „User interface“ on page 46.

Start device(s)

Do not start connected devices until the control software on the PC is started. Observe the instructions and notes from the operating instructions supplied with the individual device!

Following the prescribed power-on sequence (start control software, then turn on devices) ensures that all devices are properly detected.

Communication

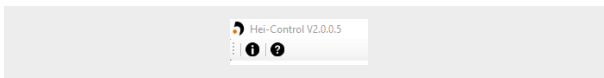
After starting and setting up the control software and switching on the connected devices, the connection is established by assigning the interfaces used via the Hei-CONTROL interface, see section „Process parameters screen“ on page 47.

User interface

After the software starts, a welcome screen appears for a few seconds, followed by the home screen.


Home screen

At the top left of the home screen menu bar, the current version of the program is displayed:


- Click the info button to open the stored manufacturer information.
- Click the question button to open the stored operating instructions in the standard browser.

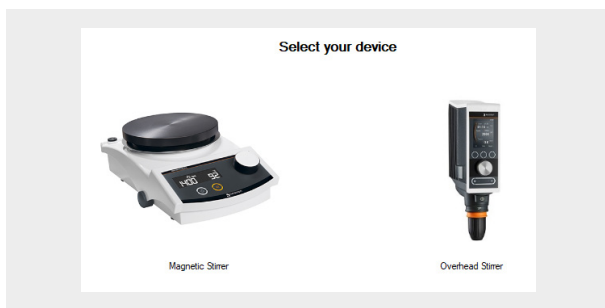The device selection appears in the main window:

- Select the desired device here by clicking on it.

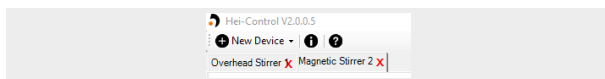In the following view, the menu bar is expanded by the **New Device** button and the tabs for the devices created (in the example, a overhead stirrer and a magnetic stirrer):

- Click the **New Device** button to add another device.
- Select the device type from the drop-down list (overhead stirrer or magnetic stirrer).
 - One tab is added per device.
 - The associated **Process Parameters** screen is opened automatically when a device is created (see section „Process parameters screen“ on page 47).
 - Each device entry can be deleted by clicking on the corresponding tab symbol [X]. Confirm the prompt to remove it.

Process parameters screen

In the **Process Parameters** screen, you can define the process parameters for a device.

The screen is opened automatically after a new device has been created, see section „Home screen“ on page 46:

The image shows two side-by-side screenshots of the 'Process Parameters' screen. The left screen is for a 'Laboratory stirrer' and the right screen is for a 'Magnetic stirrer'. Both screens have a black background with white text and input fields. The 'Laboratory stirrer' screen has sections for 'Vessel' (Size, Diameter), 'Medium' (Name, Volume, Medium temperature, Ambient temperature), 'Impeller' (Type, Size, Material), and 'Note'. The 'Magnetic stirrer' screen has sections for 'Vessel' (Size, Diameter), 'Medium' (Name, Volume, Medium temperature, Ambient temperature), 'Stirring bar' (Size, Shape), and 'Note'. Both screens have a 'Create' button and a 'Skip >' button at the bottom.

- Enter the process parameters for the selected device:
 - **Vessel:** Vessel size and diameter
 - **Medium:** Name of the medium, volume and temperature, ambient temperature
 - **Impellers:** Stirring tool data (laboratory stirrer) or
 - **Stirring bar:** Stirring Bars data (magnetic stirrer)
 - **Note:** Field for user-specific notes
- Save the data record by clicking **Create**.

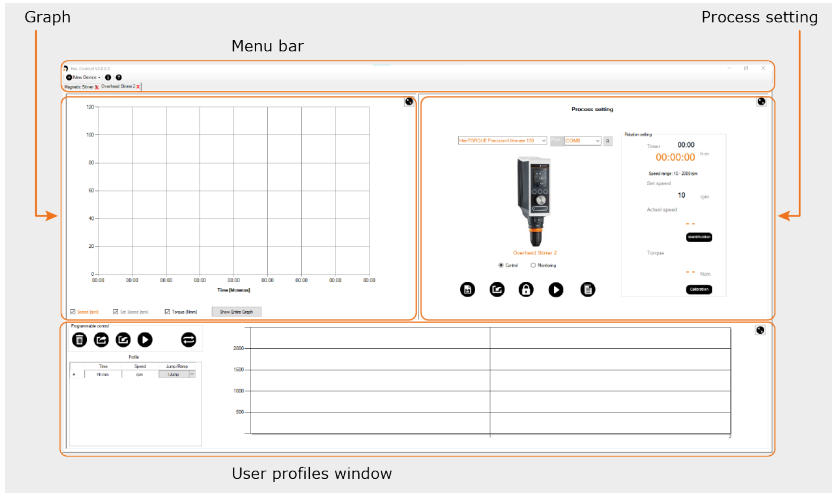
or

- Skip the mask by clicking **Skip** (the process data can be defined and edited at any time).

After confirming with **Create** or **Skip**, the process view of the last device created is opened (in the following example: overhead stirrer).

Process view

The process view is divided into three sections. The functions correspond to the menu navigation of the connected device:

The following sections describe the process view controls and input fields.

A description of the specific device functions (parameterization, handling of graphs and profiles) and parameters can be found in the operating instructions for the respective laboratory or magnetic stirrer.

Control elements

Symbol	Function
	Shortcut on the Windows® desktop to start the program.
	Zoom button: Switch between integrated view and full screen of the window area.
	Save process parameters button: Store current values as a process data file on your hard disk drive (format *.csv).
	Load process parameters button: Load stored process data file into the system.
	Save profile/process data button: Save new/modified profile or process values on your hard disk drive.
	Load profile/process data button: Load stored profile/process data.
	Lock button: Locks all controls to protect against unauthorized use.
	Unlock button: unlock all controls.
	Start button: Start loaded profile / activated device.
	Stop button: Stop the current profile / activated device.
	Delete button: Reset the loaded profile.
	Repeat button: Repeat the loaded profile.

Graphics area

Depending on the selection, the temperature and/or speed characteristics are displayed in the Graphics area.

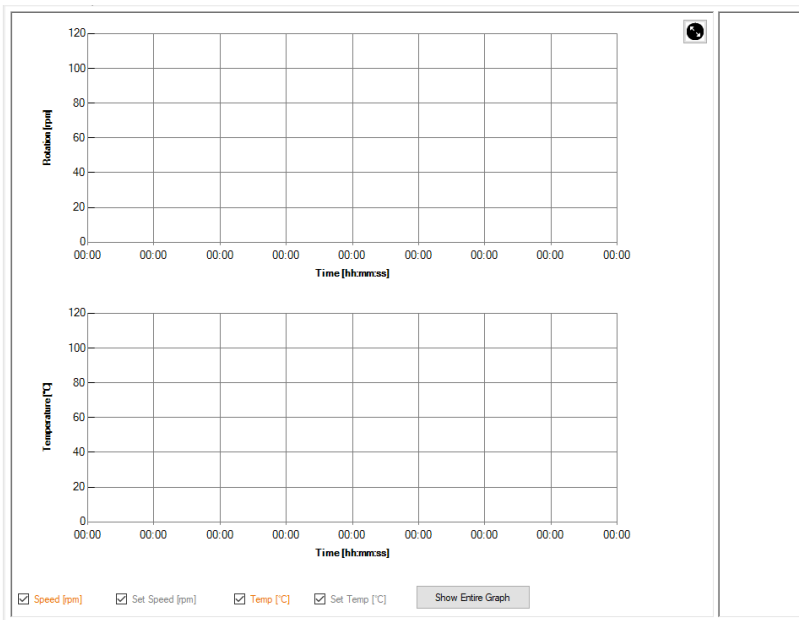
Click the zoom button in the upper right corner to switch between integrated view and full screen.

Select the check boxes at the bottom of the screen to customize the display:


- Overhead stirrer:
 - Speed [rpm]: Actual speed value
 - Set speed [rpm]: Speed setpoint
 - Torque [Nm]: Actual torque value

- Magnetic stirrer:
 - Speed [rpm]: Actual speed value
 - Set speed [rpm]: Speed setpoint
 - Temp [°C]: Actual temperature value
 - Set Temp [°C]: Temperature setpoint


Magnetic stirrer (as an example):

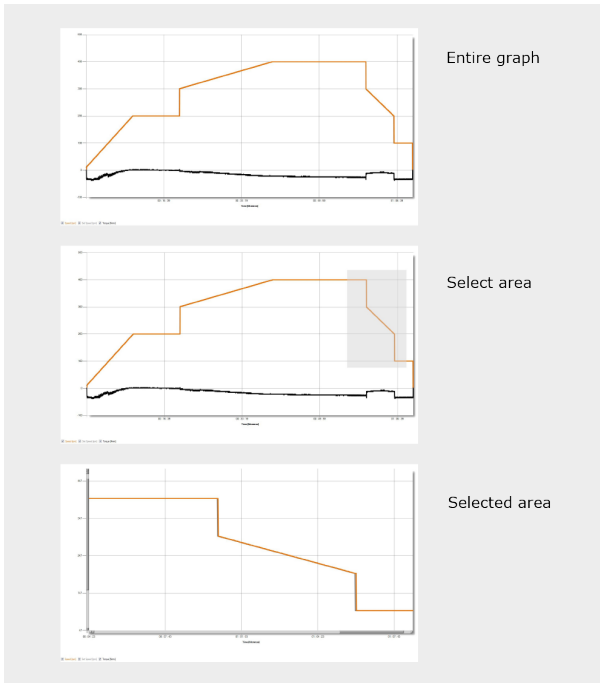Click the **Show entire Graph** button to open the Picture-in-Picture view with zoom function. This view shows the area recorded up to this point in time (example view):

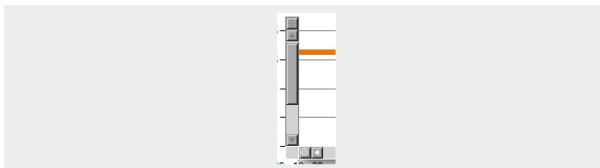
- Select a specific section (right mouse button, drag frame).
 - The selected section is zoomed in and immediately displayed.
- Click on the reset icons of the scroll bars to zoom out again:

Process parameters area

The process parameters area is used to select the device variant, assign the interface used, and select the operating mode.

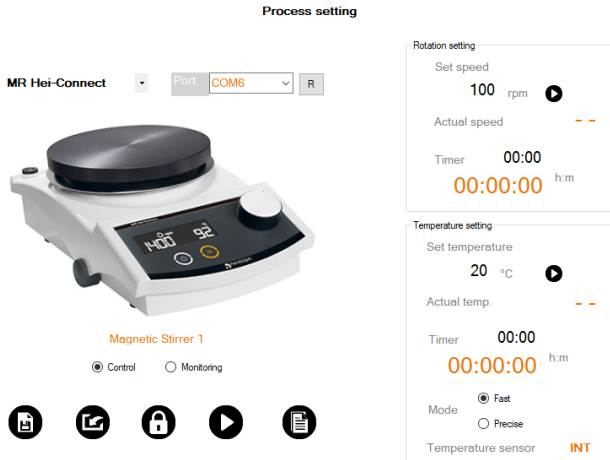In **Control** mode (see following section „Control devices“ on page 53), the process parameters speed, stirring time and temperature can be set.

The following options are available:

- Selection of the device type: In the [Device] drop-down list, select the connected device type, in the example: **MR Hei-Connect**
- Selection of the communication interface: From the [Port] drop-down list, select the interface you are using, in the example: **COM6**
- Apply device settings: Click the [R] button to accept the set values. The connection to the device can then be established.
- Load process data: Click the **Load process data** to reload stored process data.
- Lock function: Click the **Lock** button to set a password to protect against unauthorized operation/changes. Enter your valid password to unlock.

To unlock without a valid password, the tab belonging to the device must be closed, see section „Home screen“ on page 46.

- Start/stop the connected device with the **Start/Stop** buttons.
- Load process parameters: Click the **Load process parameters** button to load stored process parameters.

For a description of the controls, see section „Control elements“ on page 49.

Monitor devices

In the process parameters area, select **Monitoring** if you want to monitor connected devices using the software.

In **Monitoring** mode, all device functions are operated directly on the device.

- The parameters **Rotation setting** (magnetic and laboratory stirrers) and **Temperature setting** (magnetic stirrers only) cannot be edited. All function keys in these fields are hidden.

Control devices

In the process parameters area, select **Control** if you want to control connected devices using the software.

In **Control** mode, basic device functions are controlled by software:

- The parameters **Rotation setting** (magnetic and laboratory stirrers) and **Temperature setting** (magnetic stirrers only) can be adjusted by the user.

Value	Function
Set speed	Speed setpoint: Mark this value and enter the desired speed via the keyboard of your PC.
Actual speed	Actual speed value (read only).
Timer	Timer function: Highlight this value and enter the desired stirring time using the keyboard of your PC.
Set temperature	Temperature setpoint: Mark this value and enter the desired maximum temperature via the keyboard of your PC (magnetic stirrer only).
Actual temp.	Actual temperature value (magnetic stirrer only, read only).
Fashion	Fast / Precise heating mode selection (magnetic stirrers only, for more information on the heating modes, see the operating instructions for the activated magnetic stirrer).
Temperature sensor	Temperature sensor indicator (internal or external, magnetic stirrer only).
Torque	Actual torque value (laboratory stirrer only, read only).

- Laboratory stirrers only: The **[Range]** button can be used to select the gear stage of a connected Hei-TORQUE Precision/Ulimate 400. For device variants 100 and 200, the key designation changes to **[Identification]** and is not functional.
- Laboratory stirrers only: The **[Calibration]** button can be used to calibrate the torque. For more information on torque calibration, see the operating instructions for the controlled Hei-TORQUE.
- If desired, save the specified process parameters to reload at any time: **Save process parameters** and **Load process parameters** buttons.

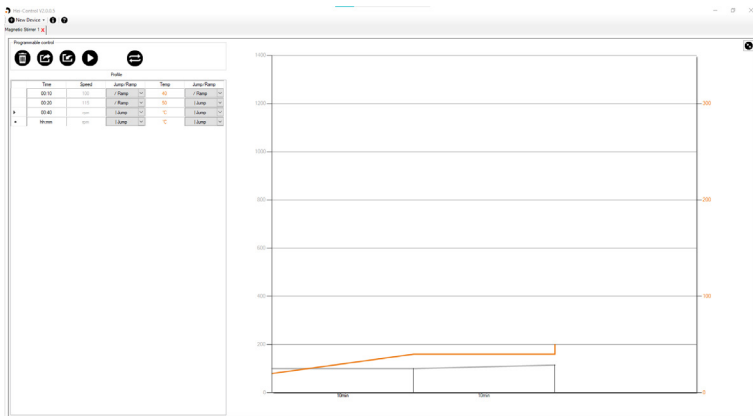
User profiles area

In the user profiles area, user-specific profiles can be created and edited in the same way as the device software:

To create a user-specific profile, proceed as described below:

- Place the cursor in the **Time** field.
 - The value is highlighted. At the same time, another input line is inserted.
- Use the keyboard to enter the stirring time value for the step.
- Place the cursor in the **Speed** field.
- Use the keyboard to enter the speed value for the step.
- In the corresponding **Jump/Ramp** drop-down list, specify whether the speed should be abruptly (jump) or gradually (ramp) augmented.
- Place the cursor in the **Temp** field.
- Use the keyboard to enter the temperature value for the step.
- In the corresponding **Jump/Ramp** drop-down list, specify whether the temperature should be abruptly (jump) or gradually (ramp) augmented.
 - Each process step that has been defined is immediately shown in the diagram:

- Define as many further process steps as required, as described.
- Start the new profile by clicking the **Start** button.
- Save the new profile by clicking the **Save profile** button.
- Load an existing profile by clicking the **Load profile** button.
- Delete a loaded profile by clicking the **Delete** button.

- Activate the repeat function by clicking the **Repeat** button.
 - The active profile is then processed repeatedly.

The cursor can be placed in the next input field with the input key and with the tab key of your PC keyboard.

The values of a loaded profile can be edited at any time.

For a description of the controls, see section „Control elements“ on page 49.

Troubleshooting

System message	Cause / possible remedy
The program has been modified or is already running!	EXE file has been renamed: Rename the .EXE file to its original name Program already started: Confirm message with OK
Connection error!	Interface cable not plugged in/loose: Connect the cable/check that it is securely attached Incorrect COM port assigned: Correct assignment
No data to save	No data entered: Enter profile information
Impossible to save profile	Save canceled: Finish saving the profile with Save
Impossible to restore profile: Wrong parameters in file	Invalid data in profile: Re-enter and save profile data
Impossible to restore profile	Undefined error loading profile: Re-enter and save profile data
Port xxx can't be opened!	Selected COM port already occupied: Assign to another COM port

Contact information Heidolph international

Heidolph Instruments North America

Phone: 1-866-650-9604
E-mail: service@heidolph.com
www.heidolphNA.com

Heidolph Instruments United Kingdom

Phone: 01799 - 5133-20
E-mail: service@radleys.co.uk
www.heidolph-instruments.co.uk

Local distributors

To find your local distributor please visit www.heidolph.com

Warranty Statement

Heidolph Instruments assumes a three-year warranty against material and manufacturing defects.

Excluded from the warranty are glass and wear parts, transport damage, and damage resulting from improper handling or non-intended use of the product.

For registered products, the warranty period begins on the date of purchase. Register the product with the enclosed warranty card or on our homepage www.heidolph.com.

For non-registered products, the warranty period begins with the date of the serial production (to be determined by the serial number).

In the event of material or manufacturing defects, the product will either be repaired or replaced free of charge within the warranty period.

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© Heidolph Instruments GmbH & Co. KG