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Zertifikate – Certifications

Contents

About this document.....	16
Typographic conventions.....	16
Copyright protection.....	16
Safety instructions in the official languages of the European Union	16
California Residents.....	16
Intended use.....	17
Compliant use.....	17
Reasonably foreseeable misuse	17
General safety instructions.....	17
Other regulations.....	17
Residual risk	17
General notes	18
Setup, Positioning	18
Magnetic stirrer – lab jack cable connection	19
Limiting the travel distance	20
Operation	21
Lowering the lab jack automatically.....	22
Signaling.....	22
Technical specifications	23
Scope of delivery	23
General cleaning instructions.....	23
Repairs - Return of equipment.....	23
Maintenance.....	24
Disposal.....	24
Registration in accordance with the Electrical and Electronic Equipment Act (ElektroG).....	24
Warranty statement.....	24
Contact details	24
Certificate of decontamination.....	25

About this document

These mounting instructions describe the operation of e-Lift type lab jacks. The mounting instructions manual is an integral part of the delivery.

Typographic conventions

This document uses the following symbols, signal words, and highlights:

	Description
	Warning symbols in combination with a signal word indicate dangers: DANGER Indicates an immediate dangerous situation. Failure to respect the indications will result in death or serious injury. WARNING Indicates a potential danger. Failure to respect the indications will result in serious injuries. CAUTION Indicates a potential hazard. If not avoided, damage to property and minor to moderate injuries can occur.
	Mandatory signs are used to indicate important and useful information on handling a product. This information is used to ensure operational safety and to maintain the value of the product.
[...]	Square brackets identify labels of control elements on the device as well as labels and entries in software masks and user interfaces.
→	The arrow indicates specific instructions to be followed to ensure operational safety when handling the product.

Copyright protection

This document is protected by copyright and is intended for use by the purchaser of the product only.

No transfer to third parties, reproduction in any form, including excerpts, and by any means, as well as utilization and/or disclosure of the contents is permitted without the prior written consent of Heidolph Scientific Products GmbH. Any violation is subject to compensation for damage.

Safety instructions in the official languages of the European Union

A summary of all safety instructions in this document, in all official languages of the European Union, can be found in our Safety Guide for the Hei-PLATE product group (Ref. 01-005-006-68). This document is available for download on our homepage in the most up-to-date version.

California Residents

Important information for California residents regarding Prop 65:
Please visit www.P65Warnings.ca.gov for more information.

Intended use

The variant of the described product delivered by Heidolph Scientific Products GmbH has been developed and designed for operation with Hei-PLATE Mix'n'Heat Ultimate magnetic stirrers (lifting and lowering of the magnetic stirrer placed on the lab jack). Any other use of this device is not considered as intended!

Compliant use

The compliance of each individual application must always be evaluated by the user. Any additional measures necessary to ensure compliance are always the responsibility of the user.

Reasonably foreseeable misuse

Additional measures may be necessary for use under conditions or for purposes deviating from the intended use, and/or specific guidelines and safety regulations must be observed. Compliance with and implementation of all relevant directives and safety measures for the respective field of application is the sole responsibility of the operator.

All risks resulting from improper use are borne solely by the operator.

The product may only be operated by authorized and instructed personnel. Training and qualification of the operating personnel as well as ensuring that the product is operated with responsibility are the sole responsibility of the operator!

General safety instructions

Before commissioning and using the product, familiarize yourself with all the safety regulations and occupational safety guidelines applicable at the place of use and observe them at all times.

Only operate the product if it is in faultless technical condition.

If there is missing or misleading information on the product or on occupational safety, contact the responsible safety specialist.

Other regulations

In addition to the notes and instructions in this document, observe all other applicable regulations such as laboratory and workplace guidelines, hazardous substances ordinances, recognized rules of safety engineering and occupational medicine as well as particular local regulations!

Noncompliance will invalidate any warranty claims against Heidolph Scientific Products GmbH.

The operating company is solely liable for all damage resulting from unauthorized changes or modifications to the device, from the use of unapproved or non-genuine spare parts and accessories, or from disregarding the safety instructions and hazard warnings or the manufacturer's instructions!

Residual risk

The product has been designed and manufactured in accordance with the state-of-the-art standards and the recognized safety regulations at the time of the development. During mounting and use, as well as during maintenance, repair and cleaning work, there are nevertheless certain residual risks associated with the described device.

These are identified and described at the appropriate points in this document.

General notes

The optional e-Lift lab jack can be used to raise and lower manually or lower automatically Hei-PLATE Mix'n'Heat Ultimate magnetic stirrers of Heidolph Scientific Products GmbH placed on the platform in laboratory operations. The lab jack can be moved not only via the integrated hand-wheel but also via the control elements and the software of the connected magnetic stirrer, see the associated operating instructions Hei-PLATE Mix'n'Heat Expert/Ultimate (01-005-006-69).

CAUTION

The USB-C interface of the magnetic stirrer may only be used for operation with a Heidolph Scientific Products GmbH lab jack. The general USB-C standard is not supported. Direct or indirect damage caused by connecting devices not authorized by the manufacturer is the responsibility of the user.


WARNING

The lab jack is designed for a load of maximum 10 kg. Direct or indirect damages caused by an overload of the lab jack, are generally the sole responsibility of the user.

Always be sure to position all setups and accessory parts centrally above the magnetic stirrer on the lab jack.

Move the lab jack downward and/or pull the setups centrally upward as much as possible to remove glass flasks or attachments from the magnetic stirrer!

Never swing setups outwards (see example: Setup with magnetic stirrer, heating attachment and Findenser®).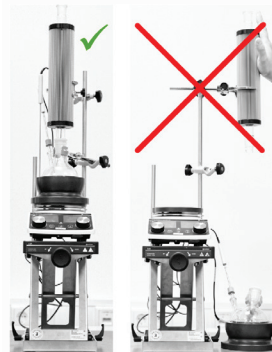

Setup, Positioning


Thanks to its robust design and the self-locking scissor mechanism, the e-Lift type lab jack supplied by Heidolph Scientific Products GmbH offers the best stability and work safety.

Follow the safety instructions below:

- To use the lab jack, place it on a clean, dry, stable, even and horizontal surface.
- Always make sure that all setup components and accessory parts of the magnetic stirrer are securely fixed (see also the corresponding delivery documentation or operating and/or mounting instructions):
- Fix the setups preferably on a wall grid (if available). When mounted on a wall grid, ensure that the wall grid and all mounting elements have sufficient self-protection and load capacity!
- Use all recommended retaining and fixing elements for setup components and accessory parts.
- The lab jack, the magnetic stirrer and any other setups do not form a physical unit and must always be transported separately.
- Before each transport, disconnect the cable connection between the lab jack and the magnetic stirrer.

Magnetic stirrer – lab jack cable connection


The lab jack [2] is connected to the USB-C connection socket at the back of the magnetic stirrer [1] using the USB-C cable supplied: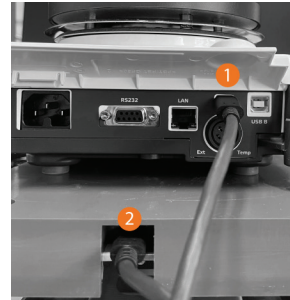

As soon as the lab jack has been detected by the software, the corresponding symbol appears on the display of the magnetic stirrer:

In case of interruption of the data connection between the lab jack and magnetic stirrer, the following warning appears on the display of the magnetic stirrer [73: Lab jack has no connection] and the [Lab jack connected] symbol goes out.

This warning message can be acknowledged at any time, and this has no negative effects on the ongoing processes of the magnetic stirrer.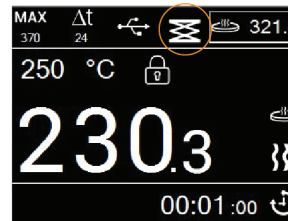

CAUTION

Please note that for ensuring the operation and process safety, the lab jack has to be permanently connected to the magnetic stirrer via the supplied USB C cable and has to be permanently supplied with power!

In the currentless state (e.g. during a power failure, during lacking cable connection between lab jack and magnetic stirrer or when the magnetic stirrer switches to standby mode), there is a risk that the lab jack lowers under high load and that processes get eventually falsified in case of an automatic restart.



If necessary, deactivate the function [Automatic restart] of the magnetic stirrer.

Limiting the travel distance


The travel distance of the lab jack can be limited mechanically with the help of suitable and adjustable clamping levers (optional accessory):

Each adjustable clamping lever [1] is mounted with a suitable shim [2] and a guide piece [3]:


- Place the guide piece [3] in the guide of the lab jack [4a] from the inside.
- Place the shim [2] on the thread of the clamping lever [1].
- Screw the clamping lever [1] onto the guide piece [3]:
 - Fix the second clamping lever in the same way in the opposite guide of the lab jack [4b].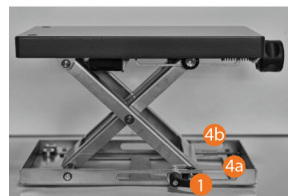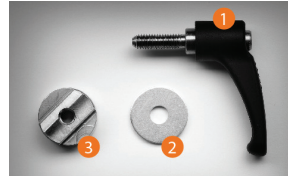

- Move the lab jack upward to the highest working point [5], push the two clamping levers onto the two guide pins [6] until they stop and tighten the two clamping levers.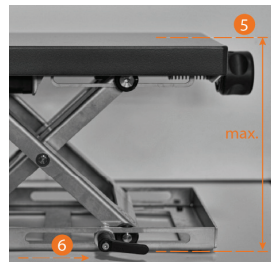

- The position of the clamping lever can be adjusted, if necessary: Pull the clamping lever outwards to unlock it [1a]. Turn the clamping lever clockwise or counterclockwise in a suitable position [1b].

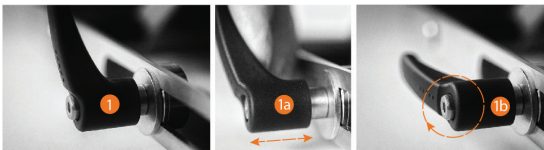The lab jack can now not be moved beyond this vertical stop point. Check both thumbscrews for tight fit before commissioning and regularly during operation.

Operation

CAUTION

- Never reach into the lab jack mechanism after it is connected to the magnetic stirrer.
- Do not place any objects in the lab jack mechanism.
- Always make sure that there is sufficient distance between the lab jack and other devices or structures: the required minimum distance is 15 cm on all sides.
- Make sure that no incoming or outgoing lines (power connection cable, data cable, hose connections) protrude into the lab jack's mechanism.
- Clean the lab jack's mechanism at regular intervals or immediately if it is highly soiled.
- Always make sure that the travel path of the lab jack is not blocked.
- In the event of high setups, pay particular attention to moving the center of gravity of the whole setup to the middle of the lab jack table as far as possible to ensure adequate stability.
- Attach a suitable warning of unexpected traversing movements of the lab jack to the overall setup in remote and in unsupervised operation.

The lab jack is moved up and down using the two control buttons (up)/(down) or the handwheel. Use the handwheel especially for fine adjustment:

- Turn the handwheel clockwise to move the lab jack downward.
- Turn the handwheel counterclockwise to move the lab jack upward.



General notes

If no lab jack is detected by the magnetic stirrer software, the information [No lift connected] appears on opening the [Lift] menu. In this case, check the cable connection between the magnetic stirrer and lab jack.

If the lab jack has been detected, the [Lift connected] symbol appears in the header of the display (first figure). As soon as the [Lower lift automatically] function is activated ([Timer expired] and/or [Error] option activated) the [Lift connected] symbol is marked with the safety symbol (second figure):

In automatic mode, the lab jack can only move downward. The upward traversing movement is only controlled via the [Lift] button and the handwheel at the front!

When lowering the lab jack, the movement speed is significantly reduced 150 mm before reaching the physically possible end position due to safety reasons.

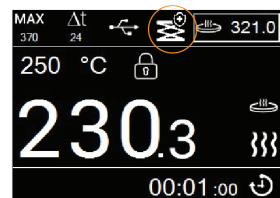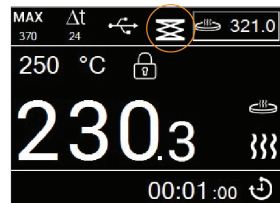
- In case of a sensor error (warning code W68), the lab jack can be moved down and up until the end position using the control buttons on the magnetic stirrer.
- As long as the sensor error exists, the downward movement speed is automatically and significantly reduced.
- If the lab jack is moved upward, the upper end position may possibly not be detected. In this case, the motor keeps turning and the click sound can still be heard until you release the button. In this case, the system cannot be damaged!
- If the sensor error occurs in automatic mode, the lab jack is moved until the lower end position with a significantly reduced speed and regardless of the set travel distance.

Lowering the lab jack automatically

From the Main menu, open the [Lift] menu and make the required settings for automatic lowering of the lab jack:

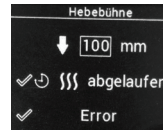- Define the required travel distance in the first line first.
- Mark the (heating) [Timer expired] option if the lab jack is to move downward automatically after the timer has expired. If the option is deselected, the lab jack does not move after the timer has expired. An eventual sensor error has no influence on the timer function.
- Mark the [Error] option if the lab jack is to move downward immediately in case of an error (error message or warning message!). The magnetic stirrer then switches to the safe operating state (speed 100 rpm, heating off). If the option is deselected, the lab jack does not move in case of error.

With automatic lowering, the lift jack is positioned with an accuracy of ± 5 mm.

Signaling

When moving the lab jack, the following visual and/or acoustic messages are generated:

- When moving the lab jack automatically, a warning symbol [Risk of crushing] is shown on the display of the magnetic stirrer for the whole moving duration.
- When moving the lab jack upward and downward, the magnetic stirrer generates a clack sound in every operating state and for the whole moving duration.
- Button sounds ON: lab jack moves automatically downward after a timer has expired: Magnetic stirrer generates two or five signal sounds.
- Button sounds OFF: lab jack moves automatically downward after a timer has expired: Magnetic stirrer generates a signal sound.
- Button sounds ON: lab jack moves automatically downward in case of an error: Magnetic stirrer generates a signal sound.
- Button sounds OFF: lab jack moves automatically downward in case of an error: Magnetic stirrer generates a signal sound.
- On reaching the lower end stop, the warning message [W71/W72] is also generated in normal operation, since this is interpreted by the software as "Motor blockage" / "End stop reached".

Technical specifications

Lab jack

Load capacity
Protection class (EN 60529)
Power supply
Footprint (W × D)
Usable top plate (W × D)
Height
Weight

e-Lift

max. 10 kg
Mechanic IP02, electronics IP42
From magnetic stirrer by USB-C cable
185 × 390 mm, incl. handwheel
185 × 350 mm
min. 95 mm, max. 275 mm
6 kg

Scope of delivery

Item	Quantity	Product no.
e-Lift lab jack	1	509-00100-00
USB type C connection cable	1	14-300-022-14
Warranty registration	1	01-006-002-78

General cleaning instructions

Wipe all surfaces and the control panel with a damp cloth if necessary. Persistent contamination can be removed with mild soapy water.

CAUTION

Clean the device's surfaces with a soft, lint-free and only slightly moistened cloth.

Never use any aggressive or abrasive cleaning agents or aids.

Repairs - Return of equipment

Repairs to the device may only be carried out by authorized experts! Unauthorized repairs during the warranty period will result in the loss of the warranty claim. Regardless of the warranty claim, the owner is solely liable for damage caused by unauthorized repairs.

- In case of repair and before returning your device, contact our technical service at the following e-mail address:
 - service@heidolph.de.
- In your message, please provide us with the following information in addition to an error description:
 - Item number
 - Serial no.

The required data can be found on the rating plate of the device.

A service representative will contact you as soon as possible to agree on the next steps.

Include the completed certificate of decontamination with every device return, see "Certificate of decontamination" on page 25.

Maintenance

The device contains no user-serviceable components. If necessary (in the event of abnormal operating behavior such as excessive noise or heat generation, for example) contact our technical service.

Disposal

- When disposing of the device, observe the provisions of the WEEE Directive 2012/19/EU and its transposition into national law in the country of use.
- When disposing of portable batteries, observe the provisions of the European Battery Directive 2013/56/EU and their implementation in national law in the country of use.
- Check the device and all components for residues of substances that are hazardous to health, the environment and biohazardous before disposal.
- Properly remove and dispose of residues of substances that are hazardous to health, the environment and biohazardous!

Registration in accordance with the Electrical and Electronic Equipment Act (ElektroG)

Heidolph Scientific Products GmbH, headquartered in 91126 Schwabach/Germany, address Walpersdorfer Str. 12, is registered under the number DE 50705753 in the electronic waste equipment register of the foundation ear (www.ear-system.de).

Warranty statement

Heidolph Scientific Products GmbH provides a three-year warranty against material and manufacturing defects.

Glass and wear parts, transportation damage, and damage resulting from improper handling or non-intended use of the product are excluded from the warranty.

The warranty period for registered products begins on the date of purchase. Register the product with the enclosed warranty card or on our homepage www.heidolph.com.

For non-registered products, the warranty period begins with the date of the serial production (to be determined by the serial number).

In the event of material or manufacturing defects, the product will either be repaired or replaced free of charge within the warranty period.

Contact details

Heidolph Scientific Products GmbH

Technical service
Walpersdorfer Str. 12
D-91126 Schwabach/Germany
Email: service@heidolph.de

Representations

To find your local Heidolph distributor please visit www.heidolph.com

Certificate of decontamination

Enclose the certificate of decontamination, duly completed, with your device return. Submissions without a certificate of decontamination cannot be processed!

CERTIFICATE OF DECONTAMINATION

IN CASE OF RETURNS

Please fill in the required fields.

Note: The sender must package the goods properly and appropriately for transport.

Heidolph Scientific Products GmbH

Walpersdorfer Straße 12
91126 Schwabach

Phone: +49 (0) 9122 9920-380

Fax: +49 (0) 9122 9920-19

E-Mail: service@heidolph.de

SENDER

Name _____ First name _____
 Company/institution _____ Department _____
 _____ Workgroup _____
 Address _____
 ZC/City _____
 Country _____ Phone _____
 Email _____

DEVICE DETAILS

Article number _____ Serial no. _____

Ticket number _____

Reason for sending in _____

Has the device been cleaned, decontaminated/disinfected? Yes No (Please mark as applicable)

If yes, which measures were carried out?

Does this device pose a risk to people and/or the environment due to the processing of substances that are hazardous to health, the environment and/or are biohazardous? Yes No (Please mark as applicable)

If yes, with which substances did the device come into contact?

LEGALLY BINDING DECLARATION

The principal/consignor is aware that they are liable to the agent/consignee for losses or damage incurred due to incomplete and incorrect information.

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Schwabach, 11. September 2023

Handwritten signature of Andreas Hahn in blue ink.

Andreas Hahn
Managing Director

Handwritten signature of Jörg Ziel in blue ink.

Jörg Ziel
Quality Manager

In accordance with UK Government guidance

UK DECLARATION OF INCORPORATION

This declaration is issued under the sole responsibility of the manufacturer,

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Model: 509-00100-00-0

Description:

Device for lifting laboratory equipment.

The object of the declaration described above is classified as Partly Completed Machinery in accordance with UK Statutory Instrument 2008 No. 1597 Supply of Machinery (Safety) Regulations 2008.

The object of the declaration described above also is in conformity with the relevant UK Statutory Instruments (and their amendments):

2008 2016 No. 1091

The Electromagnetic Compatibility Regulations 2016

2012 No. 3032

The Restriction of the Use of Certain Hazardous Substances

in Electrical and Electronic Equipment Regulations 2012

This product is supplied as an optional accessory for incorporation to a Machine Tool.

The device has to be installed in accordance with the instructions contained in the Installation Guide. The partly completed machinery must not be put into service until it has been established that the machinery into which the partly completed machinery is to be incorporated complies with the provisions of a relevant product directive as a whole. And complies with the following technical standards :

EN ISO 12100:2010, EN 61326-1:2013, EN 61010-1:2010, EN IEC 63000:2018

UK Authorised Representative (for authorities only):

ProductIP (UK) Ltd.

8. Northumberland Av. London WC2N 5BY

Signed for and on behalf of Heidolph Scientific Products GmbH
Walpersdorfer Straße 12, 91126 Schwabach / Germany

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With this declaration, we confirm (according to current knowledge) that all sold laboratory devices and components by Heidolph Scientific Products GmbH fulfill the requirements of the EU directive 2011/65/EU (RoHS) and the extension 2015/863. All devices are compatible with the requirements of the RoHS Directive for the following materials:

Max. 0,01% of weight in homogenous material for cadmium and max. 0,1% of the weight in homogeneous material for Lead, Mercury, Hexavalent chromium, Polybrominated biphenyls (PBB), Polybrominated diphenyl ethers (PBDE), Bis(2-ethylhexyl) phthalate (DEHP), Butyl benzyl phthalate (BBP), Dibutyl phthalate (DBP), Diisobutyl phthalate (DIBP).

In the case of individual assemblies, maximum concentrations maybe exceeded within the permissible exceptions of the Directive.

Schwabach, 04. September. 2024

Andreas Hahn
Managing Director

Jörg Ziel
Quality Manager

China RoHS DECLARATION OF CONFORMITY

Heidolph Scientific Products GmbH has made reasonable efforts to ensure that hazardous materials and substances may not be used in its products.

In order to determine the concentration of hazardous substances in all homogeneous materials of the subassemblies, a "Product Conformity Assessment" (PCA) procedure was performed. As defined in GB/T 26572 the "Maximum Concentration Value" limits (MCV) apply to these restricted substances:

- Lead (Pb): 0.1%
- Mercury (Hg): 0.1%
- Cadmium (Cd): 0.01%
- Hexavalent chromium (Cr(VI)): 0.1%
- Polybrominated biphenyls (PBB): 0.1%
- Polybrominated diphenyl ether (PBDE): 0.1%

Environmental Friendly Use Period (EFUP)

EFUP defines the period in years during which the hazardous substances contained in electrical and electronic products will not leak or mutate under normal operating conditions. During normal use by the user such electrical and electronic products will not result in serious environmental pollution, cause serious bodily injury or damage to the user's assets.

The Environmental Friendly Use Period for Heidolph Scientific Products GmbH products is 25 years.

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This table is created according to SJ/T 11364-2014

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Note: Table applies to all products. Some of the components or parts listed above may not be part of the enclosed product.

- O: 表示该有毒有害物质在该部件所有均质材料中的含量均在GB/T 26572规定的限量要求以下。
O: Indicates that the above mentioned hazardous substance contained in all homogeneous materials of the part is below the required limit as defined in GB/T 26572.
- X: 表示该有毒有害物质至少在该部件某一均质材料中的含量超出GB/T 26572规定的限量要求。
X: Indicates that the above mentioned hazardous substance contained in at least one of the homogeneous materials of this part is above the required limit as defined in GB/T 26572.

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Apart from the disclosures in the above table, the subassemblies are not intentionally manufactured or formulated with lead (Pb), mercury (Hg), cadmium (Cd), hexavalent chromium (CrVI), polybrominated biphenyls (PBB), and polybrominated diphenyl ethers (PBDE).

Products manufactured by Heidolph Scientific Products GmbH may enter into further devices or can be used together with other appliances .

With these products and appliances in particular, Heidolph Scientific Products GmbH will not take responsibility for the EFUP of those products and appliances.

Schwabach, 04.September.2024

Andreas Hahn
Managing Director

Jörg Ziel
Quality Manager

© Heidolph Scientific Products GmbH

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Subject to change without notice. The printed version of this document is not regularly updated. The latest issue of this document can be found by visiting our homepage.

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MATERIAL CONTENT DECLARATION FOR Heidolph Scientific Products GmbH PRODUCTS