Safety Guide

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www.heidolph.com

About this document

This document contains a summary of all safety instructions for overhead stirrers, type Hei-TORQUE in the official languages of the European Union.

This document in no way replaces the operating instructions supplied with the product! For more detailed information, please refer to the original instructions (in German) or the translation of the original instructions supplied with the product.

Typographic conventions

The following symbols, signal words and highlighting elements are used in this document:

Symbol	Signal word / explanation
	<p>Warning symbols in combination with a signal word indicate dangers:</p> <p>DANGER Indicates an immediate dangerous situation. Failure to respect the indications will result in death or serious injury.</p> <p>WARNING Indicates a potential danger. Failure to respect the indications will result in serious injuries.</p> <p>CAUTION Indicates a potential hazard which, If not avoided, damage to property and minor to moderate injuries can occur.</p>
	<p>Mandatory signs are used to indicate important and useful information on handling a product. This information is used to ensure operational safety and to maintain the value of the product.</p>

Copyright protection

This document is protected by copyright and is generally intended for use by the purchaser of the product.

No transfer to third parties, reproduction in any form, including excerpts, and by any means, as well as utilization and/or disclosure of the contents is permitted without the prior written consent of Heidolph Instruments GmbH & Co. KG. Any violation is subject to compensation for damage.

Residual risk

The device has been designed and manufactured in accordance with the state-of-the-art standards at the time of development and the recognized safety regulations. During mounting and use, as well as during maintenance, repair and cleaning work, there are nevertheless certain residual risks associated with the described device.

These are identified and described in the present document.

Intended use

The Hei-TORQUE overhead stirrers have been specifically developed for the following tasks: Stirring, mixing, gassing and degassing, emulsifying, suspending.

The Hei-TORQUE overhead stirrers are suitable for use in the following areas: Chemical, pharmaceutical, biology, environmental analysis, basic research, research laboratories.

Any other use of the described device is not considered as intended!

Due to its design, in the as-delivered condition, use in the food, cosmetics, and pharmaceutical industries as well as other comparable industries that manufacture products intended for consumption by humans or animals, or for use on humans or animals is only permitted in analytical processes or under laboratory-like conditions.

Reasonably foreseeable misuse

Additional measures may be necessary for use under conditions or for purposes deviating from the intended use, and/or specific guidelines and safety regulations must be observed. Corresponding requirements must be evaluated and implemented by the operator in each individual case.

Compliance with and implementation of all relevant directives and safety measures for the respective field of application is generally the responsibility of the operator.

All risks resulting from improper use are borne generally by the operator.

The device may generally be operated by authorized and instructed personnel. Training and qualification of the operating personnel as well as ensuring that the device is handled responsibly are the general responsibility of the operator!

Compliant use

The user is generally responsible for evaluating the conformity of his application and, if necessary, for taking additional measures.

Transportation

During transportation, avoid severe shocks and mechanical stresses that can cause damage to the device. Keep the original packaging in a dry and protected place for later use.

Storage

Always store the device in its original packaging. To protect against damage and excessive material aging, store the device in an environment that is as dry, temperature-stable and dust-free as possible.

Acclimatization

After each transportation and after storage under critical climatic conditions (e.g. high temperature difference between inside and outside), allow the device to acclimatize at room temperature for at least two hours to prevent possible damage due to condensation before commissioning it in the place of use. If necessary, extend the acclimatization phase if the temperature differences are very high.

Make all supply connections (power supply, tubing) only after the device has been acclimatized!

Permissible ambient conditions

The device is designed for indoor use only. The device is not suitable for outdoor use! The device is not suitable for use in potentially explosive areas!

When used in corrosive atmospheres, the service life of the device may be reduced depending on the concentration, duration and frequency of exposure.

General safety instructions

Before commissioning and using the device, familiarize yourself with all the safety regulations and occupational safety guidelines applicable at the place of use and observe them at all times.

Only operate the device if it is in faultless technical condition. In particular, make sure that the device and the supply connections are free of visible damage.

If there is missing or misleading information on the device or regarding occupational safety, contact the responsible safety specialist or our technical service.

Only use the device in accordance with the regulations for intended use.

Electrical safety

Before connecting the device to the power supply, ensure that the voltage indicated on the rating plate matches the specifications of the local power utility company.

Ensure that the power supply circuit provided is protected by means of a residual-current device (RCD).

Always use the power supply cord provided with the device.

Prior to use, check that the device and the power supply cord are free of visible damage.

Have repairs and/or maintenance work on the device carried out exclusively by an authorized electrician or by the technical service department of Heidolph Instruments.

Always disconnect the device from the power supply system before carrying out any maintenance, cleaning or repair work.

Data security

The user is responsible for ensuring data security when transferring data between the described device and other devices.

Always use secure networks for the data transfer and avoid use of critical infrastructure.

Always use high-quality shielded data cables for the data transfer.

Operational safety

Operate the device under a closed ventilated fume hood when working with potentially hazardous substances (see EN 14175 and DIN 12924).

Do not make any unauthorized changes or modifications to the device!

Always use genuine spare parts and accessories, or those expressly approved by the manufacturer!

Rectify malfunctions or faults on the device immediately.

Switch off and disconnect the device from the power supply, preventing reconnection, if it is not possible to eliminate the malfunction or rectify the fault immediately.

Observe all other applicable regulations such as laboratory and workplace guidelines, recognized safety technology rules and special local regulations.

Occupational safety

Always use the prescribed personal protective equipment (PPE) such as protective clothing, safety goggles, protective gloves, safety shoes, etc.

Do not operate any other devices in the immediate vicinity of the device ...

- which can generate electromagnetic fields in the frequency range between 9×10^3 Hz to 3×10^{11} Hz,
- which generate emission or radiation sources in the frequency range 3×10^{11} Hz to 3×10^{15} Hz (in the optical spectral range wavelengths from 1,000 μm to 0,1 μm),
- which generate ultrasonic or ionizing waves.

Do not operate the unit when adiabatic compression or shock waves may occur (pressure wave ignition).

Do not use substances that could release energy in an uncontrolled way and cause a pressure increase (exothermic reaction, spontaneous ignition of dusts).

Only use the stirrer tools approved by the manufacturer.

Route all cables free of kinks and outside the operating and hazardous area.

Avoid excessive pressure on the device display.

Keep the base unit dry during operation.

Ensure adequate safety distance: Do not store objects in the working and hazardous area of the device during operation.

Operate the device under a closed ventilated fume cupboard when working with potentially hazardous substances/ materials (see EN 14175 and DIN 12924).

Personal protective equipment (PPE)

The operating company must determine and provide the necessary PPE, depending on the respective application and the media and chemicals used.

The corresponding instruction of the personnel is generally within the operating company's responsibility.

Environmental protection

When processing environmentally hazardous substances, take appropriate measures to avoid hazards to the environment.

The evaluation of corresponding measures such as the marking of a hazardous area, their implementation, and the training of the relevant personnel is the general responsibility of the operator!

Biohazard

When processing biohazardous substances, take appropriate measures to prevent hazards to persons and the environment, including:

- Instruction of the personnel regarding the necessary safety measures.
- Provision of personal protective equipment (PPE) and instruction of the personnel in its use.
- Marking the device with a biohazard warning symbol.

The evaluation of corresponding measures such as the marking of a hazardous area, their implementation, and the training of the relevant personnel is the general responsibility of the operator!

Other regulations

In addition to the notes and instructions in this document, observe all other applicable regulations such as laboratory and workplace guidelines, hazardous substances ordinances, recognized rules of safety engineering and occupational medicine as well as particular local regulations!

Noncompliance will invalidate any warranty claims against Heidolph Instruments.

The operator is generally liable for all damage resulting from unauthorized changes or modifications to the device, from the use of unapproved or non-genuine spare parts and accessories, or from disregarding the safety instructions and hazard warnings or the manufacturer's instructions!

Specific safety instructions

Display

CAUTION

Pointed or sharp objects can damage the surface of the display on the control panel!

To operate the device functions, always touch the display with your fingertips or a special stylus with a soft-touch tip.

Interfaces

WARNING

If the voltages at the connection sockets are too high and insulation is inadequate, voltage can be applied to metallic parts such as the housing in case of a fault.

Extra-low voltage inputs and outputs above 25 V AC or 60 V DC respectively must be disconnected safely according to EN 61140 or must be protected by doubled or increased insulation in accordance with EN 60730-1.

Only use shielded connection cables and connect the shielding to the connector housing.

Make sure that the interface connections at the back of the device are secured by the covers provided.

Set up the device

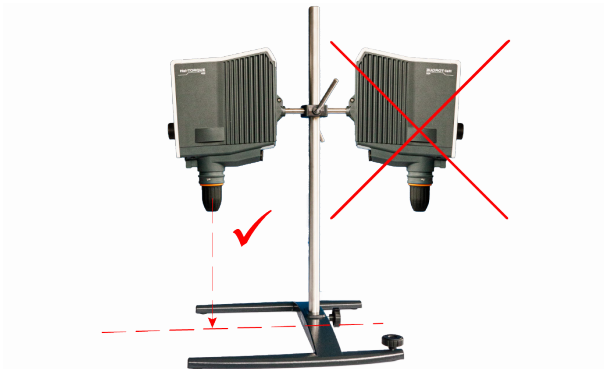
WARNING

If assembled improperly, the device can also fall or tip over when unloaded!

If the overhead stirrer is mounted on a wall grid, ensure that the wall grid and all mounting elements have sufficient self-protection and load capacity!

When mounting the overhead stirrer on a stand, make sure that the setup is sufficiently stable.

- The stand may only be set up on a stable surface.
- When mounting on a stand, follow the specific setup instructions for the stand.
- Always align the overhead stirrer so that the quick-action chuck vertically faces the notional centerline between the **long** stand base bars, see following figure:

After assembly of the overhead stirrer, check all screw connections / clamping screws for tight fit.

Quick-action chuck (Quick Chuck)

WARNING

Before mounting/dismantling the quick-action chuck, switch off the device and disconnect it from the power supply to prevent unexpected starting-up of the device!

Clamping/removing stirrer tools

WARNING

If necessary, stop the rotational movement and before clamping/removing stirrer tools, switch off the device to prevent it from starting up unexpectedly!

Automatic restart

CAUTION

Always activate or deactivate the automatic restart function as required before the actual process start!

Pay attention to the safety symbols on the device for signaling the functional status for all process sequences.

If the automatic restart function is active, secure the device with an information sign if necessary.

Starting/stopping rotation

WARNING

Risk of injury due to glass break and leaking medium

- Select the diameter of the laboratory vessel so that the stirrer tool turns inside the vessel without touching it.
- After clamping the stirrer tool and before the process starts, check the stirrer tool and the chuck for smooth running.
- Never start up stirrer tools or chucks that have an unbalance, and replace them immediately.

Risk of injury on rotating parts

- Never grip turning parts.
- Always wear close-fitting clothing when working with the device.
- Do not wear any accessories such as ties, scarves, shawls, pieces of jewelry, etc. when working with the device.
- When working with the device, tie back long hair or wear a cap or hair net.

Damage to property, loss of production

- Adjust the start-up behavior of the device to the specific properties of the sample to be processed to avoid overflowing/ejection of the sample or overloading of the stirrer tool (safety settings).
- Follow the specific safety instructions for the substances to be processed (refer to the safety data sheets if applicable).
- Always wear the required individual personal protective equipment (PPE).

Remote mode

WARNING

Secure the device in remote mode with a clearly visible warning sign and, if necessary, take further ambient-specific protective measures that protect against damage to property and injuries in the event of unexpected/unmonitored starting of the device.

Device service

DANGER

Switch the device's main switch off and disconnect it from the power supply before carrying out maintenance work, cleaning, or repairs.

When cleaning, avoid the penetration of liquids.

Before replacing the fuses, switch off the device and disconnect the power supply.

Always replace the two device fuses in pairs with original manufacturer fuses. Further information on available accessories can be found on our website at www.heidolph.com! After fuse replacement, check the device for a safe condition according to IEC 61010-1.

CAUTION

Clean the device's surfaces with a soft, lint-free and only slightly moistened cloth.

Never use any aggressive or abrasive cleaning agents or aids.

Repairs

Repairs to the device may always be carried out by authorized experts! Unauthorized repairs during the warranty period will result in the loss of the warranty claim. The owner is generally liable for damage caused by unauthorized repairs. If repairs are necessary, contact an authorized dealer or our technical service.

Maintenance

There are no user-serviceable components in the unit housing. If necessary (in the event of abnormal operating behavior such as excessive noise or heat generation, for example) contact our technical service.

Check the caps and fasteners of all glass components for correct and tight fit at regular intervals and especially before each start-up of the device.

Check the level sensor in the heating bath regularly for proper operation, contamination, and limescale build-up to prevent malfunctions.

Disposal

When disposing of the device, observe the provisions of the WEEE Directive 2012/19/EU and its transposition into national law in the country of use.

When disposing of portable batteries, observe the provisions of the European Battery Directive 2013/56/EU and their implementation in national law in the country of use.

Check the device and all components for residues of substances that are hazardous to health, the environment and biohazardous before disposal.

Properly remove and dispose of residues of substances that are hazardous to health, the environment and biohazardous!

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© Heidolph Instruments GmbH & Co. KG