Betriebsanleitung
Operating instructions
Notice d'instructions

Hei-SHAKE Orbital Expert

 **heidolph**
research made easy

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Seite 4 – 37

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Schnittstellenbefehle RS232 (ASCII)

Generic commands

Send	Receive	Description
"IDENT\x\n"	"< Name > < MaterialNr > < SerialNr > < Fertigungsauftrag > \x\n"	Read identification numbers
"SW_VERS\x\n"	"MC X1.Y1.Z1:X2.Y2.Z2, HMI A1.B1.C1:A2. B2.C2\x\n"	Read all software versions
"STATUS\x\n"	"STATUS X\x\n"	Read actual status
"IN_SP_ALL\x\n"	"SP_3 = X, SP_10 = Y, SP_11 = Z\x\n"	Read all setpoints
"IN_PV_ALL\x\n"	"PV_5 = X, PV_12 = Y S\x\n"	Read all actual values
"CC_ON\x\n"	"CC_ON\x\n"	Set connection control
"CC_OFF\x\n"	"CC_OFF\x\n"	Reset connection control
"RESET\x\n"	"RESET X\x\n"	Reset communication

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Translation of the original instructions
Page 42 – 75

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Contents

Introduction

About this document.....	42
Typographic conventions.....	42
Copyright protection.....	42
Safety instructions in the official languages of the European Union	42

General product information

Directives applied, product certification.....	43
Copyright notice	43
Residual risk	43
California Residents.....	43
Intended use.....	43
Compliant use.....	44
Reasonably foreseeable misuse	44
Transportation.....	44
Storage.....	44
Acclimatization	44
Permissible ambient conditions	44
Energy saving potential and cost efficiency	44

Safety

General safety instructions.....	45
Electrical safety.....	45
Interfaces.....	45
Data security	45
Operational safety	46
Occupational safety	46
Electrostatic charging of liquid or powdered media.....	46
Personal protective equipment (PPE).....	46
Environmental protection.....	47
Biohazard.....	47
Other regulations.....	47

Device description

Mechanical design– control panel– terminals.....	48
--	----

Commissioning

Positioning the device	49
Power supply	49
Automatic restart	50
Setup	51
Starting/deceleration ramp	51
Reset to factory settings	51
Device parameters	52
Decimal separator	52
Display brightness.....	52
Date, Time	52
Display lock	52
Error list	52
Info.....	52
Placing and attaching laboratory vessels	53
Safety functions	53
Speed limitation	53
Display lock	54
Device behavior when the display lock is activated.....	54
Start inhibit	54

Operation

Switching the device on/off	55
Operation via touch display	55
Remote control status indicator	56
Home screen	56
Setting the shaking frequency	57
Timer function	58
Operating modes.....	59
Continuous operation	60
Timer mode.....	60
User profiles	61
Create/modify user profiles	61
Processing user profiles	62
Delete user profiles	63
Remote mode.....	64
Data logging.....	65

Troubleshooting

Troubleshooting.....	67
Warning codes	68
Error codes	68

Appendix

Technical specifications	69
Performance range/maximum load	70
RS232 (ASCII) interface commands	71
Scope of delivery	72
Accessories	72
Device service.....	72
General cleaning instructions.....	72
Repairs - Return of equipment.....	73
Maintenance.....	73
Disposal.....	73
Registration in accordance with the Electrical and Electronic Equipment Act (ElektroG)	73
Warranty statement.....	74
Contact details	74
Certificate of decontamination.....	75

About this document

This operating manual describes the features and operation of platform shakers Hei-SHAKE Orbital Expert type.

The operating instructions are an integral part of the described device!

Typographic conventions

This document uses the following symbols, signal words, and highlights:

	Description
	Warning symbols in combination with a signal word indicate dangers: DANGER Indicates an immediate dangerous situation. Failure to respect the indications will result in death or serious injury. WARNING Indicates a potential danger. Failure to respect the indications will result in serious injuries. CAUTION Indicates a potential hazard. If not avoided, damage to property and minor to moderate injuries can occur.
	Mandatory signs are used to indicate important and useful information on handling a product. This information is used to ensure operational safety and to maintain the value of the product.
[...]	Square brackets identify labels of control elements on the device as well as labels and entries in software masks and user interfaces.
→	The arrow indicates specific instructions to be followed to ensure operational safety when handling the product.

Copyright protection

This document is protected by copyright and is intended for use by the purchaser of the product only.

No transfer to third parties, reproduction in any form, including excerpts, and by any means, as well as utilization and/or disclosure of the contents is permitted without the prior written consent of Heidolph Scientific Products GmbH. Any violation is subject to compensation for damage.

Safety instructions in the official languages of the European Union

A summary of all safety instructions in the official languages of the European Union can be found in our Safety Guide for the product group Hei-SHAKE (shakers and mixers, Ref. 01-005-006-99). This document is available for download on our homepage in the most up-to-date version.

Directives applied, product certification

CE marking

The device meets all requirements of the following directives:

- Machinery Directive 2006/42/EC
- EMC Directive 2004/108/EC
- RoHS 2 directive 2011/65/EU, including amendment 2015/863

Copyright notice

The software implemented in this product is protected by copyright laws. The rights holder is Heidolph Scientific Products GmbH, Walpersdorfer Straße 12, 91126 Schwabach/Germany. Any open source software components in this software are excluded from our copyright. Further information is available in the service area on our website www.heidolph.com.

Violations of copyright (e.g. unauthorized use or modification of the software) may result in civil claims (e.g. omission, compensation for damages) and/or entail criminal penalties within the meaning of the legislation of the destination country.

Residual risk

The device has been designed and manufactured in accordance with the state-of-the-art standards at the time of development and the recognized safety regulations. During mounting and use, as well as during maintenance, repair and cleaning work, there are nevertheless certain residual risks associated with the described device.

These are identified and described at the appropriate points in this document.

California Residents

Important information for California residents regarding Prop 65. Please visit www.P65Warnings.ca.gov for more information.

Intended use

The described platform shaker in this document has been specially developed for the following tasks:

- Shaking
- Mixing
- Emulsifying
- Suspending
- Separating
- Dissolving
- Staining

The areas of application of the described platform shaker include chemical, biological and environmental-analytical laboratory and research applications, applications of basic research and comparable facilities.

Due to its design, the device in its delivery condition may only be used in analytical processes or in laboratory-like conditions in the food, cosmetics, and pharmaceutical industries as well as other comparable industries that manufacture products intended for consumption by humans or animals, or for use on humans or animals.

Any other use of this device is not considered as intended!

Compliant use

The compliance of each individual application must always be evaluated by the user. Any additional measures necessary to ensure compliance are always the responsibility of the user.

Reasonably foreseeable misuse

Additional measures may be necessary for use under conditions or for purposes deviating from the intended use, and/or specific guidelines and safety regulations must be observed. Corresponding requirements must be evaluated and implemented by the operator in each individual case.

Compliance with and implementation of all relevant directives and safety measures for the respective field of application is the sole responsibility of the operator.

All risks resulting from improper use are borne solely by the operator.

The device may only be operated by authorized and instructed personnel. Training and qualification of the operating personnel as well as ensuring that the device is handled responsibly are the sole responsibility of the operator!

Transportation

During transportation, avoid severe shocks and mechanical stresses that can cause damage to the device. Keep the original packaging in a dry and protected place for later use.

Storage

Always store the product in its original packaging. To protect against damage and excessive material aging, store the device in an environment that is as dry, temperature-stable and dust-free as possible.

Acclimatization

After each transportation and after storage under critical climatic conditions (e.g., high temperature difference between inside and outside), allow the device to acclimatize at room temperature for at least two hours to prevent possible damage due to condensation before commissioning it in the place of use. If necessary, extend the acclimatization phase if the temperature differences are very high.

Make all supply connections only after the device has been acclimatized!

Permissible ambient conditions

The device is designed for indoor use only. The device is **NOT** suitable for outdoor use! The device is **NOT** suitable for use in potentially explosive areas!

When used in corrosive atmospheres, the service life of the device may be reduced depending on the concentration, duration and frequency of exposure.

Energy saving potential and cost efficiency

Even in standby mode, small amounts of energy are used to supply individual device functions. However, especially over longer periods of nonuse, this resource consumption adds up and causes avoidable costs.

Therefore, always switch off the device when not in use, especially for long periods of time (devices with main switch: Main switch in position [O]; devices without main switch: Unplug) to effectively avoid unnecessary energy consumption and the resulting costs.

General safety instructions

Before commissioning and using the device, familiarize yourself with all the safety regulations and occupational safety guidelines applicable at the place of use and observe them at all times.

Only operate the device if it is in faultless technical condition. In particular, ensure that there is no visible damage on the device itself and, where applicable, on connected devices or the supply connections.

If there is missing or misleading information on the device or regarding occupational safety, contact the responsible safety specialist or our technical service.

Only use the device in accordance with the regulations for intended use.

Electrical safety

Before connecting the device to the power supply, ensure that the voltage indicated on the rating plate matches the specifications of the local power utility company.

Ensure that the power supply circuit provided is protected by means of a residual-current device (RCD).

Always use the power supply cord provided with the device.

Prior to use, check that the device and the power supply cord are free of visible damage.

Make sure that the power plug can be reached directly at any time, in order to disconnect the device from the power supply at any time without delay.

Have repairs and/or maintenance work on the device carried out exclusively by an authorized electrician or by the technical service department of Heidolph Scientific Products GmbH.

Always disconnect the device from the power supply system before carrying out any maintenance, cleaning or repair work.

Interfaces

Safely separate extra low voltage inputs and outputs rated 25 V AC or 60 V DC or above according to EN 61140, or use double or reinforced insulation according to EN 60730-1 or DIN 60950-1.

Use only shielded connection cables. Connect the shield to the connector housing. Always cover unused interfaces (front and back) to prevent liquids from entering the device.

Data security

The user is responsible for ensuring data security when transferring data between the described device and other devices.

Only use secure networks for the data transfer and avoid use of critical infrastructure.

Only use high-quality shielded data cables for the data transfer.

For data transfer via USB, an industrial standard USB hub should be preferably used to ensure the most stable connection possible.

Operational safety

Operate the device under a closed ventilated fume hood when working with potentially hazardous substances (see EN 14175 and DIN 12924).

Do not make any unauthorized changes or modifications to the device!

Only use genuine spare parts and accessories, or those expressly approved by the manufacturer!

Rectify malfunctions or faults on the device immediately.

Switch off and disconnect the device from the power supply, preventing reconnection, if it is not possible to eliminate the malfunction or rectify the fault immediately.

Observe all other applicable regulations such as laboratory and workplace guidelines, recognized safety technology rules and special local regulations.

Occupational safety

Always use the prescribed personal protective equipment (PPE) such as protective clothing, safety goggles, protective gloves, safety shoes, etc.

Do not operate any other devices in the immediate vicinity of the device ...

- which can generate electromagnetic fields in the frequency range between 9×10^3 Hz to 3×10^{11} Hz,
- which generate emission or radiation sources in the frequency range 3×10^{11} Hz to 3×10^{15} Hz (in the optical spectral range wavelengths from 1,000 μm to 0.1 μm),
- which generate ultrasonic or ionizing waves.

Do not process any substances that could release energy in an uncontrolled manner (e.g. self-ignition).

Do not process substances in which the energy input through mixing poses a danger.

Wipe off any liquid that may have spilled on the device immediately.

Before each transfer or transport (e.g., removal from the fume hood, transfer between laboratory benches, temporary storage in the laboratory area, etc.), always remove all attachments from the shaking platform of the device.

Electrostatic charging of liquid or powdered media

Liquid or powdered media as well as the laboratory vessels used can become dangerously electrostatically charged due to the shaking movement or when emptying or cleaning the vessel. The intensity of the charge generated depends on the properties of the medium, the flow rate, the process parameters as well as the size, geometry and material of the laboratory vessels (source: TRGS 727, Section 4).

The evaluation of a potential explosion hazard due to electrostatic charging and the definition of suitable safety measures are always the responsibility of the operator and are always based on the substances to be processed and the individual application.

Personal protective equipment (PPE)

The operating company must determine and provide the necessary PPE, depending on the respective application and the media and chemicals used.

The required instruction of the personnel is solely within the operating company's responsibility.

Environmental protection

When processing environmentally hazardous substances, take appropriate measures to avoid hazards to the environment.

The evaluation of appropriate measures such as the marking of a hazardous area, their implementation, and the training of the relevant personnel is the sole responsibility of the operator!

Biohazard

When processing biohazardous substances, take appropriate measures to prevent hazards to persons and the environment, including:

- Instruction of the personnel regarding the necessary safety measures.
- Provision of personal protective equipment (PPE) and instruction of the personnel in its use.
- Marking of the device with the biohazard warning symbol.

The evaluation of appropriate measures such as the marking of a hazardous area, their implementation, and the training of the relevant personnel is the sole responsibility of the operator!

Other regulations

In addition to the notes and instructions in this document, observe all other applicable regulations such as laboratory and workplace guidelines, hazardous substances ordinances, recognized rules of safety engineering and occupational medicine as well as particular local regulations!

Noncompliance will invalidate any warranty claims against Heidolph Scientific Products GmbH.

The operating company is solely liable for all damage resulting from unauthorized changes or modifications to the device, from the use of unapproved or non-genuine spare parts and accessories, or from disregarding the safety instructions and hazard warnings or the manufacturer's instructions!

Mechanical design– control panel– terminals

- 1 Shaking platform
- 2 Touch panel with display
- 3 [Standby] button

RS232 interface (Sub-D, 9-pin)

- 4 The RS232 interface of the device can be used for maintenance and service purposes (e.g. firmware update) as well as for controlling the device and for reading out device data. Observe the instructions in section "Remote-Betrieb" on page 17 and "RS232 (ASCII) interface commands" on page 71.
- 5 USB-A interface
- 6 USB-B interface
- 7 Appliance inlet: Power supply connection
- 8 Rating plate: Device and manufacturer data, serial number (example).

Positioning the device

CAUTION

Improper positioning of the device may result in direct and indirect damage to property due to falling and/or spilling of fluids!

The surface must be even, non-slip, proper and sufficiently stable.

The device must be freely accessible at all times.

Make sure that a sufficient safety distance to moving components can be maintained on all sides.

Observe the generally valid rules for the safe positioning of work equipment.

The professional mounting and positioning of the device including the supplied accessories is within the sole responsibility of the operator!

Heidolph Scientific Products GmbH shall under no circumstances be liable for any personal injury or damage to property resulting from failure to observe the instructions for the correct installation and positioning of the device.

Power supply

DANGER

Observe the instructions in section "General safety instructions" on page 45, in particular Electrical safety.

Power supply cord with locking system

Only use the supplied power supply cord with locking system!

Connect the power supply coupling cord [2] to the IEC- appliance inlet on the back of the device [1]. Make sure that the locking system engages [3] correctly.

Connect the power supply cord to a properly secured power socket.

Disconnecting the power supply cord

Disconnect the power supply cord from the mains socket-outlet.

Press in the locking lever on the cable coupling and disconnect the power supply cord from the IEC appliance plug on the back of the device.

Automatic restart


In the as-delivered condition, the [Automatic restart] function is disabled and has to be enabled manually.

Activate the device from standby mode. Tap the [Automatic restart] button on the Home screen menu bar, and then tap the orange icon in the selection box. Touch the button again to close the selection box.

When the [Automatic restart] function is activated (the icon in the menu bar changes to orange), the device switches on automatically upon return of the power supply after a power interruption and the speed is augmented to the last set value.

If the function remains deactivated, the device must be activated by the operator from standby mode when the mains voltage returns.

To deactivate the function, proceed in the same way: Tap the white icon in the selection field. Automatic restart is deactivated if the icon in the menu bar is displayed in white.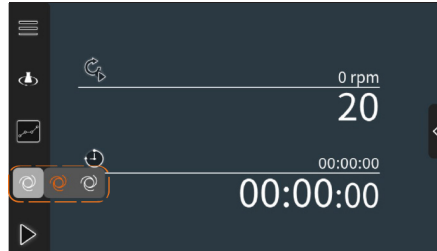

WARNING


Before activating the function, take appropriate measures to make sure that, in case of a possible unmonitored restart, no hazard for persons or risk of damage to property can occur.

The evaluation of potential risks by a possible unmonitored automatic restart as well as of appropriate security measures is the sole responsibility of the user!

Operating behavior when the [Automatic restart] function is activated

When the voltage returns after a mains voltage interruption, a warning signal sounds and a warning symbol appears on the display for five seconds. After this warning time has expired, the speed is increased again to the previously set value! The timer is restarted after each mains power interruption.

The automatic restart can be aborted within the warning time (button [Cancel] below the warning message).

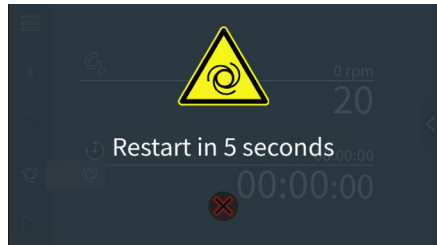Setup

Touch the [Setup] button on the Home screen menu bar to access the following settings and functions (the layout follows the menu structure of the device):

[Setup] button

[Home] button – This button takes you back to the home screen from any view.

Menu [Start/deceleration ramp settings, device reset]:

Starting/deceleration ramp

This parameter defines the starting and deceleration behavior of the device when the shaking movement is switched on or off. Options:

- [Fast]: The shaking frequency is increased or decreased as quickly as possible to the set value.
- [Medium]: The shaking frequency is increased or decreased at normal speed to the set value.
- [Slow]: The speed is gradually increased or decreased to the set value.
- [60.00 rpm/sec]: Tap this value to open the on-screen keyboard and enter a individual value. The shaking frequency is increased or decreased by the defined rate to the set value.

The setting remains active even if the power supply is interrupted!


In general, the actual start-up behavior depends on other parameters such as the total load and the load distribution on the shaking platform. The ideal setting must be determined and defined by the user in individual cases.

Reset to factory settings

Touch the [Reset] button to restore the factory settings.

Confirm the security prompt with [Yes] to permanently reset all device and operating parameters, or cancel the operation with [No].

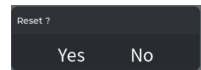Device parameters

Decimal separator

Use this selection to specify the separator for the representation of decimal numbers.
Options: Point, comma


Display brightness

Use the slider to set the desired brightness.

Date, Time

Tap the button to display the current setting: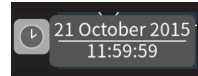

Tap the [Date] field to open the input screen for setting the system date:

Tap the [Time] field to open the input screen for setting the system time:

Display lock

Tap the button to activate the display lock. The icon changes from white to orange. In addition, the [Lock Display] icon is permanently displayed on the home screen.

To disable the display lock, tap the button again. The icon returns to white, and the [Lock Display] icon disappears from the Home screen.

See section "Device behavior when the display lock is activated" on page 54 for detailed information.

Error list

Tap the button to display the list of all recorded errors, see section "Troubleshooting" on page 67 for more information.

Info

Touch this button to open the info menu with the following options: Serial number, HMI version, micro-controller version, QR code for further information and technical documents.

Placing and attaching laboratory vessels

For placing and attaching your laboratory vessels, use the various Heidolph attachments which are available as an option. For detailed information on the available accessories, see www.heidolph.com → Products → Platform shakers → Accessories, as well as the Heidolph Assembly Instructions Hei-SHAKE, Ref. 01-005-007-13.

At low shaking frequencies, vessels such as Petri dishes can also be placed without using any kind of attachment on the non-slip contoured rubber mat.


- Always place individual vessels in the center of the shaking platform.
- Always distribute several vessels evenly on the shaking platform.

Safety functions

Speed limitation

Touch the [Speed Limit] button to open the selection to set a maximum and/or minimum shaking frequency or an RCF value.

Tap the desired value to open the on-screen keyboard. The input is made using the keys or the slider, see previous section "Setting the shaking frequency" on page 57.

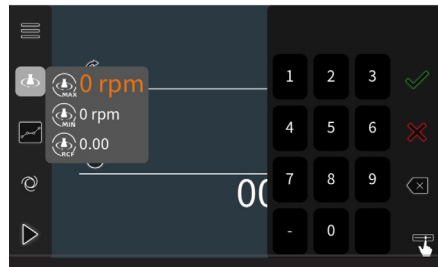If a maximum and/or minimum speed is specified, values above or below the defined limit value can no longer be entered.

When a relative centrifugal acceleration (RCF) value is set, the instrument automatically determines a corresponding speed. The maximum adjustable RCF value is 6.2 and must be selected according to the speed limits of the device.

Unsaved entries are discarded without a safety warning.

If no input is made for about twelve seconds in edit mode or if the input value is not accepted within this time, the display jumps back and the start screen is displayed again.

Display lock

To prevent unwanted operation by unauthorized persons or by accidental touching or droplets on the touch screen, it is recommended that the display lock be activated in correspondingly hazardous areas or in the case of appropriate application (e.g. sample processing with open containers).

In the as-delivered condition, the display lock is deactivated .
Open the [Setup] → [Device Parameters] menu and tap the [Display Lock] button to activate the function.

The button icon changes from white to orange, the display is locked, the touch function is deactivated:

Exception: The [STANDBY] button can be used to put the device into standby mode even if the display is locked!

When the display lock is active, the [Lock Display] icon (1) is permanently displayed on the home screen.



Device behavior when the display lock is activated

- If the lock function is active, the display is automatically locked if no operation is performed for a period of 100 seconds.
- Lock the display immediately at any time by tapping the [Lock Display] icon (1) on the Home screen.

In both cases, the lock icon and the unlock slider appear on the display.


Drag the slider to the right to unlock the display. No further entries (e.g. password) are required.

To disable the display lock, navigate to the [Device Parameters] menu again and tap the [Display Lock] button. The button icon returns to white and the [Lock Display] icon disappears.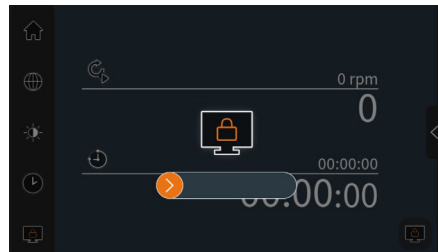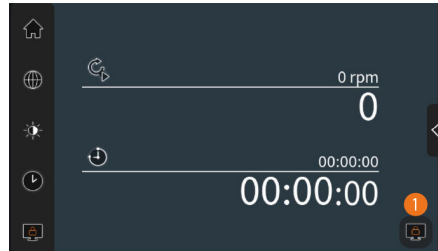

Start inhibit

To prevent an uncontrolled start-up, a start-up inhibit has been implemented in the software of the device: After starting the shaking movement with one of the [Start] or [Start profile] buttons (see following section "Home screen" on page 56), the operation must be released by dragging the slider (see figure on the right).

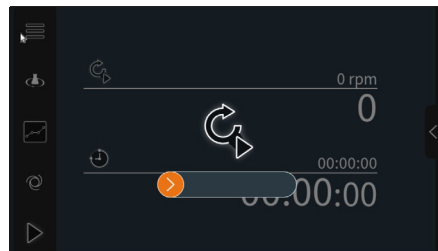Switching the device on/off

When the mains voltage is applied, the device is in standby mode (the [STANDBY] button LED lights up white). Briefly press the [STANDBY] button to activate the device: The button LED changes to permanent green.

The manufacturer logo appears briefly on the display, followed by the start screen.

Then press the [STANDBY] button to briefly put the device into standby mode. "Shutting down" is briefly displayed on the display, and the LED of the button changes back to white.

After pressing the [STANDBY] button, the device is booted or switched off within two to three seconds.

Unplug the device to de-energize it! When de-energized, the [STANDBY] button LED goes out.

Operation via touch display

The device features a touch display which is used to control the operational functions. Follow the safety instructions below:

CAUTION

The touch-sensitive display of the device may also react in the event of contact with liquids (drip formation, splash water, etc.).

Furthermore, droplets, fluid accumulation or condensation can restrict the touch function of the display.

This can result in faults, uncontrolled device behavior, unexpected operation and/or undesirable parameter changes.

- Always keep the display free of dirt, fluid buildup and/or condensation. Clean the glass surface of the display preferably when the display lock is activated, otherwise there is a risk of unintentionally adjusting the settings. For further cleaning instructions, see also section "Device service" on page 72
- Always activate the display lock during operation (see section "Display lock" on page 54).
- Confirm entries via the slider immediately to activate the display lock again (see section "Setting the speed - "Adjusting values without the keyboard" on page 57).

Remote control status indicator

When the device is connected to a remote controller via the RS 232 interface, the remote status is indicated on the home screen as follows:

- Remote icon gray: device is controlled locally, no input signal.
- Remote icon magenta-coloured: device operates in remote mode and is controlled by input signals.


Home screen

The following display and controls are available on the home screen:

Menu bar with buttons

- (1a) – Setup (see section "Setup" on page 51)
- (1b) – Speed limitation (see section "Speed limitation" on page 53)
- (1c) – User Profiles (see section "User profiles" on page 61)
- (1d) – Automatic restart (see section "Automatic restart" on page 50)
- (1e) – Start/Stop – button for starting and stopping the shaking movement (additional release via slider, see section "Start inhibit" on page 54).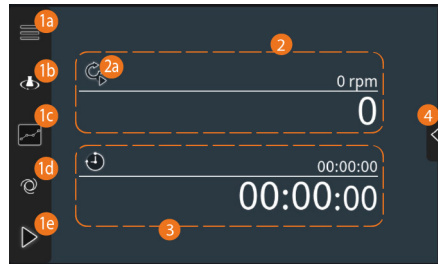

Function button Speed [2] with key [Profile Start] (2a); for more information, see sections "Start inhibit" on page 54 (additional release via slider) and "Setting the shaking frequency" on page 57.

Function button Timer [3].

Slider to hide and show the timer function button [4]. See section "Timer function" on page 58 for detailed information.

Setting the shaking frequency

In the value display area (2a) of the Speed function button (2), tap the display.

The on-screen keyboard (a) appears.

Adjustment range -530 to 530 rpm

In the negative range (-530 to -1 rpm), the shaking platform rotates clockwise.

In the positive range (1 to 530 rpm), the shaking platform rotates counter-clockwise.

The setting range can be restricted by the safety function [Speed limitation], see Section "Speed limitation" on page 53.

Function keys

To accept entries, tap [Confirm] (b).

Tap [Discard] (c) to close the on-screen keyboard without accepting the last input.

Tap [Delete] (d) to delete the last digit you entered (correct the entry).

Tap the [Slider] icon (e) to display the following options:




Adjusting values without the keyboard

Touch and drag the slider (e) up or down to adjust the speed continuously.

Touch the [+] / [-] (f) buttons to increase or decrease the speed in increments of ten.

Tap [0] (g) to set the speed to zero.

Tap the [Keyboard] icon (h) to redisplay the on-screen keyboard.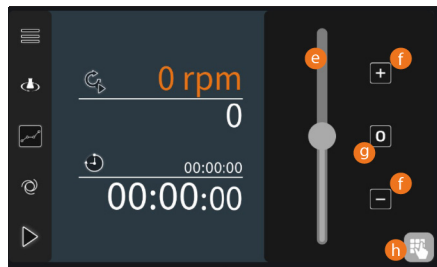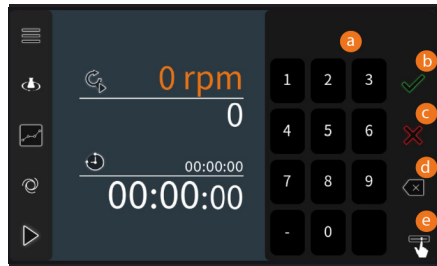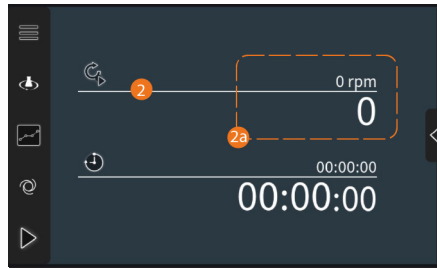

Unsaved entries are discarded without a safety warning.

If no input is made for about twelve seconds in edit mode or if the input value is not accepted within this time, the display jumps back and the start screen is displayed again.

Displaying the speed values

The upper value corresponds to the setting value in [rpm]. In edit mode, the value is displayed in large font and in orange digits. In normal operation in small font and white.

The lower value corresponds to the actual speed in [rpm]. This value is continuously updated during operation, and the display (white digits, large font) remains unchanged.

Timer function

Disabling the timer

In normal condition, the function button is displayed (timer activated, see figure on the right). Tap the slider [4] to display the [Timer] icon (orange = Timer enabled).

Tap the [Timer] icon in the slider to disable the function (icon changes to white, timer is disabled).

Tap the slider again to hide the [Timer] icon. Alternatively, the slider closes automatically after approximately three seconds.

Follow the same procedure to reactivate the timer.

Timer setting

In the value display area (3a) of the timer function button (3), tap the display.




The on-screen keyboard (a) appears.

The value [hours] is highlighted in orange: Use the on-screen keyboard to enter the desired value.

Input options

Continuous input: After three digits (e.g. 123 hours) have been entered, the display with the next digit changes to the following setting value [minutes], from this value to [seconds] after another two digits have been entered.

In this mode, leading zeros must be entered, for example: Desired runtime: 23 hours, 6 minutes and 10 seconds; input: 0230610.

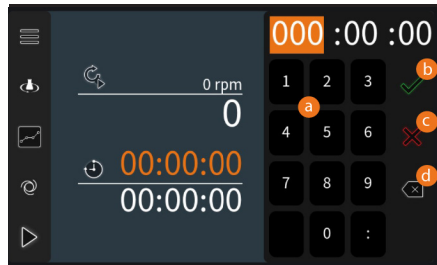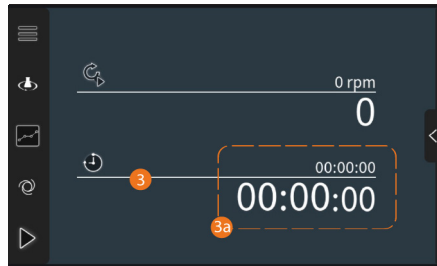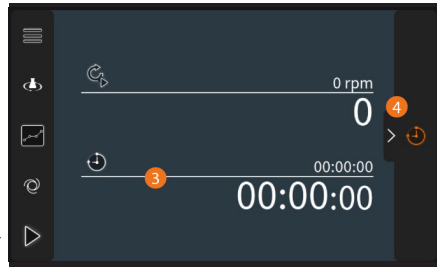Field-by-field input: Tap the desired input field [hours], [minutes] or [seconds] and enter the desired value with or without a leading zero(s).

Tap the next desired input field or use ":" to switch to the following field.

Adjustment range

The setting range of the timer function is between one second (display 000:00:01) and up to 999 hours 59 minutes and 59 seconds.

Function keys

Tap [Confirm] (b) to accept the entry.

Tap [Discard] (c) to close the on-screen keyboard without accepting the last input.

Tap [Delete] (d) to delete the last digit you entered (correct the entry).

Displaying the timer values

The upper value corresponds to the setting value in [hh:mm:ss] or [hhh:mm:ss]. In edit mode, the value is displayed in large font and in orange digits. In normal operation in small font and white.

The lower value corresponds to the time remaining until shaking is stopped in [hh:mm:ss] or [hhh:mm:ss]. This value is continuously tracked during operation, and the display remains unchanged.

The selected setting [Hide/Show Timer Function Button]. Remains active until the operator switches over again, even in the event of a power interruption.

Unsaved entries are discarded without a safety warning.

If no input is made for about twelve seconds in edit mode or if the input value is not accepted within this time, the display jumps back and the start screen is displayed again.

The (programmed) timer is started on starting the shaking movement. The function cannot be interrupted! The shaking movement can be stopped at any time before the timer expires. In this case, the last timer value remains displayed and is reset only when the shaking movement is restarted (restart timer).

Operating modes

The device can be used in continuous operation, timer controlled or remote controlled. Regardless of the selected operating mode, observe the following safety instructions:

WARNING

- Whenever possible, use closed vessels for processing corrosive, toxic or biohazardous substances and seal them safely.
- Increase the speed only gradually to the desired shaking intensity, especially with open vessels, and observe the fluid movements.
- Lower the shaking intensity or reduce the total load on the platform when the device starts to vibrate during operation.
- Always ensure that the vessels are evenly distributed on the platform.
- Pay particular attention to an even distribution of the weight load when vessels of different sizes and/or differently filled vessels are placed on the platform at the same time!
- Before switching on the device, make sure that all vessels are properly fixed on the platform.
- Use suitable attachments for the used vessels. Further information on available accessories can be found on our website at www.heidolph.com.
- Always use the necessary and appropriate personal protective equipment!


Continuous operation

In continuous operation, the shaking movement is started and stopped manually. Define the appropriate device parameters (see section "Setup" on page 51) and the desired speed (see section "Setting the shaking frequency" on page 57).

Tap one of the [Start] or [Profile Start] buttons (1) to start the shaking mode. Both buttons change from white to orange and to [Stop] or [Profile Stop] (2).

To stop shaking, tap either [Stop] or [Profile Stop] button.

If the [Timer] function button is displayed, the elapsed time since the start of the process is displayed here. If the maximum run time of 999 hours, 59 minutes and 59 seconds is exceeded, the value [>999] is displayed permanently until the process is switched off.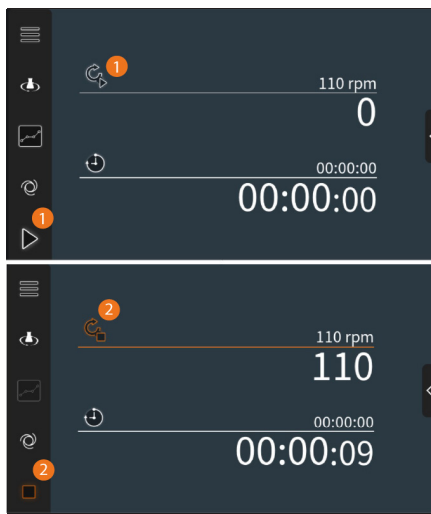

Timer mode

In timer mode, a specific period of time can be set for a continuous shaking movement.


Define the appropriate device parameters (see section "Setup" on page 51), the desired speed (see section "Setting the shaking frequency" on page 57) and the desired switch "Timer function" on page 58-on time (see section).

Tap one of the [Start] or [Profile Start] buttons (1) to start the shaking mode with timer. Both buttons change from white to orange and to [Stop] or [Profile Stop] (2). The timer icon of the [Timer] function button changes from white to orange.

If necessary, tap either [Stop] or [Profile Stop] button to stop shaking.

The remaining time until the expiration of the programmed time period is shown as [hh:mm:ss].

When the shaking mode stops, the timer is paused and will be reset when starting the shaking movement again.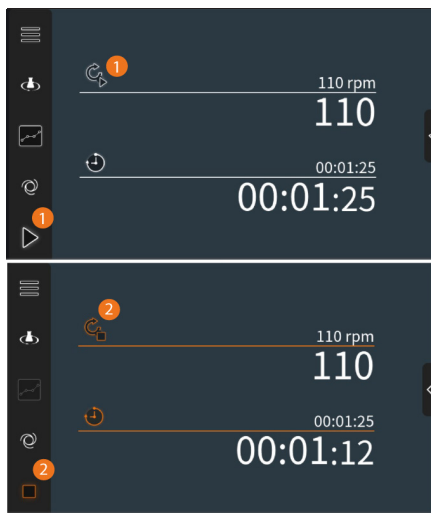


User profiles

Three user profiles, each with up to ten process steps, can be stored in the device memory.

Open the profile view using the [Profiles] key in the menu bar of the home screen (see section "Home screen" on page 56):

- (1) – Profile slots 1 / 2 / 3
- (2) – Process step indicator
- (3) – Buttons [+] and [-] to add or delete individual process steps
- (4) – Buttons [Up] and [Down] to move individual process steps
- (5) – [Delete] button to remove individual process steps

The icons of free slots are shown in white. The icons of occupied slots are displayed in orange. Used slots can be overwritten at any time.

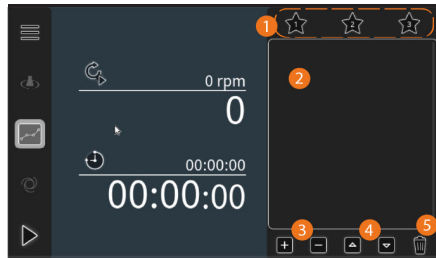Create/modify user profiles


Tap the [+] button to create a new process step (6) (in the example on the right, slot 1 - slots can be allocated as required in the last step when creating a profile).

In the [Speed] value display area, tap the display. The on-screen keyboard appears. Enter the desired shaking frequency as described in section "Setting the shaking frequency" on page 57

Tap the display in the [Timer] value display area. The on-screen keyboard appears. Enter the desired process time as described in section "Timer function" on page 58

The input values are displayed immediately in the active process step (in the example: 150 rpm for a duration of two minutes and 15 seconds):

Speed changes within a profile can occur abruptly [7] or gradually [7a].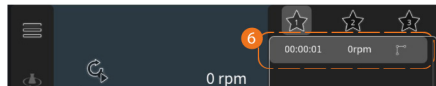

- Touch the [7]/[7a] icon of the active line to switch between [Jump] and [Ramp]: In [Jump] mode, the speed is increased or decreased as quickly as possible to the slave value. In [Ramp] mode, the speed is gradually increased or decreased to the following value.
- Create up to ten process steps per profile in the manner described.
- Move existing process steps as desired using the [Up] and [Down] arrow keys.
- Use [-] to remove unnecessary or faulty process steps.


Operation

Assign new or modified profiles to the desired slot:

Touch the icon of the desired slot for about one second until the [Save] and [Delete] icons appear (8).

Briefly tap [Save] to save the entries as a profile or to apply changes to an existing profile.

Tap [Delete] briefly to remove a profile that has already been saved.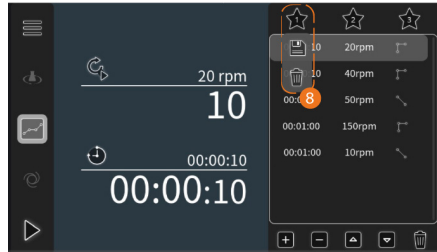

Unsaved changes are discarded when changing to another profile and exiting the view without a security prompt!


Processing user profiles

Open the profile view, briefly touch the icon of the desired slot to load the user profile.

Touch the [Start] button (9) to process the profile from process step 1.

While the profile is being processed, the [Stop] key (10) is displayed instead of the [Start] key (9).

With [Stop] or [Profile Stop], the processing of the profile can be interrupted at any time.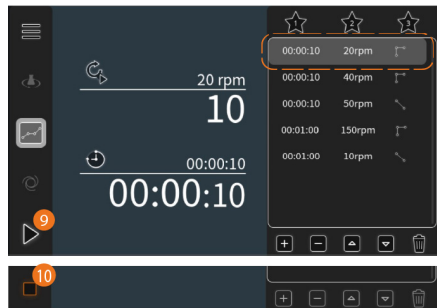

User profiles can also be partially processed, i.e. from a certain process step:

Load a profile as described above.


Select the process step from which the profile is to be processed (in the example on the right: Step 4, 150 rpm for 60 minutes).

In this case, start processing the profile with the [Profile Start] button (11).

While the profile is being processed, the corresponding stop buttons [Stop] and [Profile Stop] are displayed instead of the two start buttons.

The processing of the profile can be interrupted at any time with both stop buttons.

To cancel the process, touch the [Stop] button and exit the Profile view.

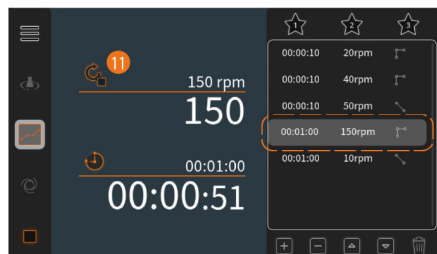Delete user profiles

Delete an individual user profile

Open the profile view, touch the icon of the assigned slot for about one second until the [Save] and [Delete] icons appear below the icon. Tap [Delete] to remove the stored profile.

Delete all user profiles

Perform a device reset. All user data is deleted, see section "Reset to factory settings" on page 51.

Remote mode

WARNING

Secure the device in remote mode with a clearly visible warning sign and, if necessary, take further ambient-specific protective measures that protect against damage to property and injuries in the event of unexpected/unmonitored starting of the device.

To activate the remote access of the device, a message of type "SetTargetSpeed" or "SetMotorState" has to be sent. Afterwards, the remote symbol appears on the display and remote commands are executed normally.

The transmission of the messages is made via ASCII strings. The protocol is active after the start. By a defined ADIP "Connect" command, you can switch to the ADIP mode.

RS232 interface parameters

- 9600 Baud
- Parity: none
- Data: 8 Bit
- Stop: 1 Bit

Command termination

All ASCII strings have to be finished with `\r\n` (0x0D 0x0A):

- ...`\r\n`

For further information on interface handling, please contact our technical service, see "Contact details" on page 74.

Device activation via PC

The device can be addressed via the integrated RS232 interface via an external control. The data interface enables activation of the device and recording of the process data in remote mode. A prerequisite for this is a suitable terminal program or software solution such as the Hei-PROCESSING Solutions from Heidolph Scientific Products GmbH

Device activation via server-based software

The device can be addressed via the integrated RS232 interface via a server-based software. If required, contact an authorized dealer or our technical service, see section "Contact details" on page 74.

Operation behavior in remote operation

In remote mode, the individual menus of the device control can be opened as described and the parameters/values can be read out, but not changed.

It is also not possible to increase or decrease the rotational speed intensity on the device (push-and-rotary control without function).


If necessary, remote operation can be terminated by the operator by pressing the [Standby] button.

Data logging

The Datalogging function is available as soon as a USB stick on the USB-A interface of the device is detected and the [USB] (1) icon is displayed on the display.

At the same time, the [Start] button is marked with a recording character (period, 2), as shown in the figure on the right.

The Datalogging function can be activated in all operating modes.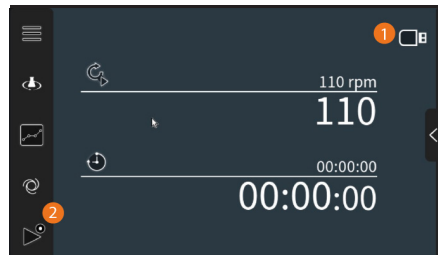

Setting the recording period



In the menu [Setup] - [Settings start/deceleration ramp, device reset], the menu item [Datalogging] (3) is also displayed for setting the desired recording duration.

Tap the [Datalogging] menu item to open the selection window for setting the recording interval (time between two data points) (4).

Options:

- off or
- 1 – 18 h or
- 1 – 59 minutes or
- 1 – 59 s

Tap the [Datalogging] menu item again to close the selection window.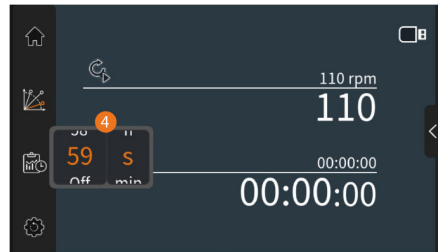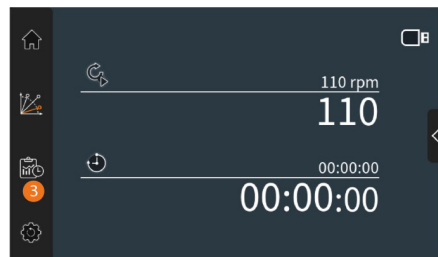

The icon of the [Datalogging] menu item is displayed in orange (5) as soon as a recording duration has been set (function activated).


When you select [off], the function remains deactivated and the icon remains white.

Start/stop datalogging


The data recording starts with the start of the shaking motion: [Start] button with white recording character switches to [Stop] button with red recording character.

Regardless of the specified recording duration, data recording is stopped if the shaking movement is interrupted by one of the two [Stop] or [Profile Stop] buttons (6).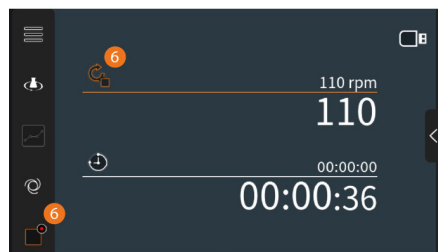

The recorded data is automatically written to the connected USB stick. During the writing process, the [USB] icon is displayed in orange.

Do not remove the USB key from the device until the icon returns to white. Otherwise, data loss may occur.

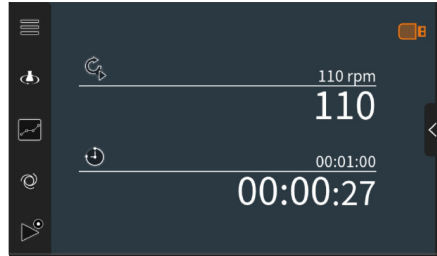Notes on how to save the device data to a USB stick

Always use USB sticks from well-known manufacturers such as Platinum TWS, XLYNE Swing, Kingston DataTraveler®, Transcend JetFlash® etc. to ensure the necessary data security.

Format the USB stick with FAT32. Use the built-in Windows function of your PC or special software.

Other formatting and USB sticks with a memory capacity greater than 32 GB are not supported in the current version of the device.

Troubleshooting

Failure	Cause	Possible remedy
Display remains dark after switching on	No mains voltage	Check the power supply cord for damages/correct connection
	Power supply cord defective	Check building fuses
	Display defective	Contact Heidolph Sales or Service
The [STANDBY] button LED does not light up when the voltage is present	Button LED defective	Contact Heidolph Sales or Service
Shaking movement stops	Motor overheat protection has tripped	Wait about 20 minutes and reduce the load on the shaking platform
	Electrical fault (no engine noise)	Contact Heidolph Sales or Service
	Mechanical defect (engine noise audible)	Contact Heidolph Sales or Service
Values "jump" during operation, speed increases or decreases unexpectedly	Strong drop formation on the touch screen	Activate the display lock
The device does not respond or is delayed to user input via the touch screen		Dry the display, if necessary Reposition superstructures on the shaking platform

In case of recurring errors please contact the responsible sales department or our technical service (see section "Contact details" on page 74).

Warning codes

Code	Description	Device behavior, remedy
W100B	Permissible PCB temperature exceeded	The device changes to a safe operating state - let it cool down, restart the process.
W100A	Permissible motor temperature exceeded	
W100C	Overvoltage	Device changes to a safe operating state - restart the process.
W100D		
W100E		
W100F		
W1010	Undervoltage	Device changes to a safe operating state - restart the process.
W1011		
W1012		
W1013		
W1014		
W1015	Engine blockage	Device changes to a safe operating state - restart the process.

Error codes

Code	Description	Device behavior, remedy
E8001	Internal memory error	Device changes to safe operating state - restart the device.
E8006	Motor error	
E8007		
E8008		
E8009	Undervoltage	Unit switches off, contact technical service.
----	HMI communication error	


- Pending error messages are displayed in orange font.
- Acknowledged error messages are displayed in white font.

Technical specifications

Hei-SHAKE Orbital Expert

Dimensions (W × H × D)	356 × 129 × 403 mm
Weight	Approx. 14 kg with perforated platform 1000 Approx. 16.5 kg with perforated platform 2000 (optional)
Interface(s)	1 × RS232 (Sub-D, 9 pole) 1 × USB A 1 × USB B
Permissible load	Max. 15 kg with perforated platform 1000 Max. 12.5 kg with perforated platform 2000 (optional)
Usable area (W × D)	294 × 263 mm with perforated platform 1000 390 × 340 mm with perforated platform 2000 (optional)
Motion	orbital vibrating
Orbit	20 mm
Speed range	-530 – 530 rpm
Drive	EC motor
Protection class	IP32 (acc. to EN 60529)
Protection class	I Ⓢ
Acoustic pressure	< 85 dB (A), in accordance with IEC 61010
Rated voltage	1/N/PE 100-240V AC 50/60 Hz
Power input	90 W
Operating and storage temperature	5 °C – 31 °C at up to 80 % rel. humidity 32 °C – 40 °C at up to 50 % rel. humidity (decreasing linearly)
Installation altitude	up to 2,000 m asl

Performance range/maximum load

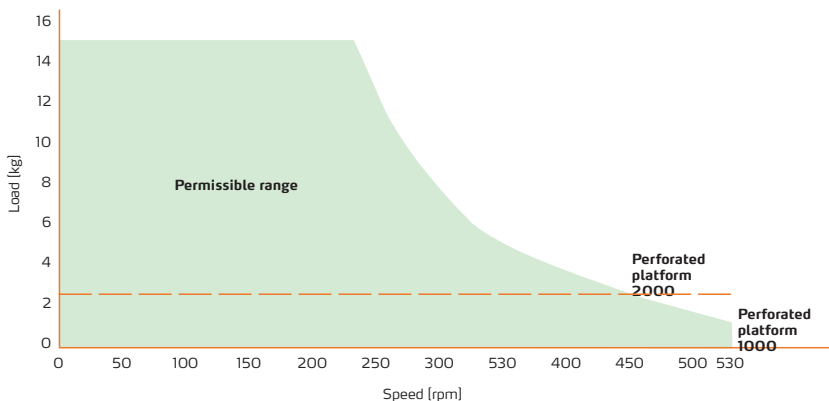The displayed values were determined with flat weight blocks and without further superstructures on the shaking platform (perforated platform 1000) and are only used for orientation.

Under these test conditions, the maximum speed when using the standard perforated platform 1000 is 530 rpm, decreasing with increasing load. When using an optional perforated platform 2000, the maximum speed is approximately 450 rpm due to the higher weight of the platform, decreasing with increasing load.

The actual maximum possible speed can be determined by the user in individual cases depending on the process and the individual design (perforated platform 1000 or 2000, brackets used, number of vessels, sample quality). It is not possible to give a general manufacturer recommendation for the different scenarios due to the application-specific parameters.

RS232 (ASCII) interface commands

Generic commands

Send	Receive	Description
"IDENT\x\n"	"< Name > < MaterialNr > < SerialNr > < Production Order > \x\n"	Read identification numbers
"SW_VERS\x\n"	"MC X1.Y1.Z1:X2.Y2.Z2, HMI A1.B1.C1:A2. B2.C2\x\n"	Read all software versions
"STATUS\x\n"	"STATUS X\x\n"	Read actual status
"IN_SP_ALL\x\n"	"SP_3 = X, SP_10 = Y, SP_11 = Z\x\n"	Read all setpoints
"IN_PV_ALL\x\n"	"PV_5 = X, PV_12 = Y S\x\n"	Read all actual values
"CC_ON\x\n"	"CC_ON\x\n"	Set connection control
"CC_OFF\x\n"	"CC_OFF\x\n"	Reset connection control
"RESET\x\n"	"RESET X\x\n"	Reset communication

Speed control commands

Send	Receive	Description
"OUT_SP_3 X\x\n"	"OUT_SP_3 X\x\n"	Set target speed (rpm)
"OUT_SP_12 X\x\n"	"OUT_SP_12 X\x\n"	Set maximum motor moment
"IN_PV_5\x\n"	"IN_PV_5 X\x\n"	Get actual speed (rpm)
"OUT_SP_14 0\x\n"	"OUT_SP_14 0\x\n"	Set Minimum Speed in rpm
"OUT_SP_15 100\x\n"	"OUT_SP_15 100\x\n"	Set Maximum Speed in rpm
"START_2\x\n"	"START_2\x\n"	Start movement
"STOP_2\x\n"	"STOP_2\x\n"	Stop movement
"OUT_MODE_10 0\x\n"	"OUT_MODE_10 0\x\n"	Select speed controller mode

Position control commands

Send	Receive	Description
"OUT_SP_10 X\x\n"	"OUT_SP_10 X\x\n"	Set target position (degrees)
"OUT_SP_12 X\x\n"	"OUT_SP_12 X\x\n"	Set maximum motor moment
"OUT_SP_13 X\x\n"	"OUT_SP_13 X\x\n"	Set maximum speed for positioning (rpm)
"IN_PV_12\x\n"	"IN_PV_12 X\x\n"	Get actual position / Get absolute position
"START_2\x\n"	"START_2\x\n"	Start movement
"STOP_2\x\n"	"STOP_2\x\n"	Stop movement
"OUT_MODE_10 1\x\n"	"OUT_MODE_10 1\x\n"	Select position controller mode

Scope of delivery

Item	Quantity	Product no.
Hei-SHAKE Orbital Expert	1	546-31400-00
Power supply cord	1	country specific
Warranty registration	1	01-006-003-16

Accessories

Further information on available accessories can be found on our website at www.heidolph.com.

Device service

When carrying out service work on the device (cleaning, maintenance, repair), observe the general instructions and safety information described in this section.

DANGER

Switch the device's main switch off and disconnect it from the power supply before carrying out maintenance work, cleaning, or repairs.

When cleaning, avoid the penetration of liquids.

Before replacing the fuses, switch off the device and disconnect the power supply cord.

Always replace the two device fuses in pairs with original manufacturer fuses. Further information on available accessories can be found on our website at www.heidolph.com.

After fuse replacement, check the device for a safe condition according to IEC 61010-1.

General cleaning instructions

Wipe all surfaces and the control panel with a damp cloth if necessary. Persistent contamination can be removed with mild soapy water.

CAUTION

Clean the device's surfaces with a soft, lint-free and only slightly moistened cloth.

Never use any aggressive or abrasive cleaning agents or aids.

Repairs - Return of equipment

Repairs to the device may only be carried out by authorized experts! Unauthorized repairs during the warranty period will result in the loss of the warranty claim. Regardless of the warranty claim, the owner is solely liable for damage caused by unauthorized repairs.

- In case of repair and before returning your device, contact our technical service at the following e-mail address:
 - service@heidolph.de.
- In your message, please provide us with the following information in addition to an error description:
 - Item number
 - Serial no.

The required data can be found on the rating plate of the device.

A service representative will contact you as soon as possible to agree on the next steps.

Include the completed certificate of decontamination with every device return, see "Certificate of decontamination" on page 75.

Maintenance

The device contains no user-serviceable components. If necessary (in the event of abnormal operating behavior such as excessive noise or heat generation, for example) contact our technical service.

Disposal

- When disposing of the device, observe the provisions of the WEEE Directive 2012/19/EU and its transposition into national law in the country of use.
- When disposing of portable batteries, observe the provisions of the European Battery Directive 2013/56/EU and their implementation in national law in the country of use.
- Check the device and all components for residues of substances that are hazardous to health, the environment and biohazardous before disposal.
- Properly remove and dispose of residues of substances that are hazardous to health, the environment and biohazardous!

Registration in accordance with the Electrical and Electronic Equipment Act (ElektroG)

Heidolph Scientific Products GmbH, headquartered in 91126 Schwabach/Germany, address Walpersdorfer Str. 12, is registered under the number DE 50705753 in the electronic waste equipment register of the foundation ear (www.ear-system.de).

Warranty statement

Heidolph Scientific Products GmbH provides a three-year warranty against material and manufacturing defects.

Glass and wear parts, transportation damage, and damage resulting from improper handling or non-intended use of the product are excluded from the warranty.

The warranty period for registered products begins on the date of purchase. Register the product with the enclosed warranty card or on our homepage www.heidolph.com.

For non-registered products, the warranty period begins with the date of the serial production (to be determined by the serial number).

In the event of material or manufacturing defects, the product will either be repaired or replaced free of charge within the warranty period.

Contact details

Heidolph Scientific Products GmbH

Technical service

Walpersdorfer Str. 12

D-91126 Schwabach/Germany

Email: service@heidolph.de

Representations

To find your local Heidolph distributor please visit www.heidolph.com

Certificate of decontamination

Enclose the certificate of decontamination, duly completed, with your device return. Submissions without a certificate of decontamination cannot be processed!

CERTIFICATE OF DECONTAMINATION

IN CASE OF RETURNS

Please fill in the required fields.

Note: The sender must package the goods properly and appropriately for transport.

Heidolph Scientific Products GmbH

Walpersdorfer Straße 12
91126 Schwabach

Phone: +49 (0) 9122 9920-380

Fax: +49 (0) 9122 9920-19

E-Mail: service@heidolph.de

SENDER

Name _____ First name _____
 Company/institution _____ Department _____
 _____ Workgroup _____
 Address _____
 ZC/City _____
 Country _____ Phone _____
 Email _____

DEVICE DETAILS

Article number _____ Serial no. _____

Ticket number _____

Reason for sending in _____

Has the device been cleaned, decontaminated/disinfected? Yes No (Please mark as applicable)

If yes, which measures were carried out?

Does this device pose a risk to people and/or the environment due to the processing of substances that are hazardous to health, the environment and/or are biohazardous? Yes No (Please mark as applicable)

If yes, with which substances did the device come into contact?

LEGALLY BINDING DECLARATION

The principal/consignor is aware that they are liable to the agent/consignee for losses or damage incurred due to incomplete and incorrect information.

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hereby declare, that the product (from serial number on 200066855) designated below is in compliance with the basic requirements of all applicable EU-directives stated below with regard to design, type of model sold and manufactured by us. This certificate will be invalid if the product is modified without the prior written consent and agreement of the manufacturer.

Reax 2	541-21001-XX		
Reax top	541-10000-XX		
Reax control	541-11000-XX		
Reax 20/4	541-20004-XX		
Reax 20/8	541-20008-XX		
Reax 20/12	541-20012-XX		
Titramax 100	544-11100-XX, 544-11200-XX		
Duomax 1030	543-32105-XX, 543-32205-XX, 543-32210-XX		
Polymax 1040	543-42105-XX, 543-42110-XX, 543-42205-XX, 543-42210-XX		
Polymax 2040	542-40005-XX, 542-40010-XX		
Multi Reax	545-10000-XX		
Vibramax 100	544-21200-XX	Titramax 101	544-11300-XX
Rotamax 120	544-41200-XX	Vibramax 110	544-31200-XX
Titramax 1000	544-12200-XX	Inkubator 1000	549-90010-XX
Unimax 1010	543-12310-XX	Promax 1020	543-22332-XX
Unimax 2010	542-10020-XX	Promax 2020	542-20020-XX
Hei SHAKE Orbital Core	546-11XXX-XX		
Hei SHAKE Salto Core	546-17XXX-XX		
Hei SHAKE ROCK	546-134XX-XX, 546-154XX-XX		
Hei SHAKE WAVE	546-144XX-XX, 546-164XX-XX		
Hei SHAKE REZIPROK	546-124xx-xx		
Hei SHAKE Orbital Expert	546-314xx-xx		

Maschinenrichtlinie / Machinery Directive 2006/42/EG
EMV-Richtlinie / Electromagnetic Compatibility Directive 2014/30/EU
Delegierte (RoHS-) Richtlinie / Delegated (RoHS) Directive 2015/863/EU
Angewandte (harmonisierte) Normen / (Harmonized) Standards applied:
EN ISO 12100:2010, EN 61326-1:2013, EN 61010-1:2010, EN IEC 63000:2018

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Andreas Hahn
Managing Director

Andreas Hahn
Managing Director

Jörg Ziel
Director Quality Management
and Technical Service

Jörg Ziel
Director Quality Management
and Technical Service

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**UK
CA**

Declaration of Conformity

In accordance with UK Government guidance

This declaration is issued under the sole responsibility of the manufacturer,

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Description:

Shaker & Mixer with different working directions

The object of the declaration described above is in conformity with the relevant UK Statutory Instruments (and their amendments):

2008 No. 1597

The Supply of Machinery (Safety) Regulations

2008 2016 No. 1091

The Electromagnetic Compatibility Regulations 2016

2012 No. 3032

The Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic Equipment Regulations 2012

and complies with the following technical standards :

EN ISO 12100:2010, EN 61326-1:2013, EN 61010-1:2010, EN IEC 63000:2018

UK Authorised Representative (for authorities only)

ProductIP (UK) Ltd.

8, Northumberland Av.

London WC2N 5BY

Signed for and on behalf of Heidolph Scientific Products GmbH

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With this declaration, we confirm (according to current knowledge) that all sold laboratory devices and components by Heidolph Scientific Products GmbH fulfill the requirements of the EU directive 2011/65/EU (RoHS) and the extension 2015/863. All devices are compatible with the requirements of the RoHS Directive for the following materials:

Max. 0,01% of weight in homogenous material for cadmium and max. 0,1% of the weight in homogeneous material for Lead, Mercury, Hexavalent chromium, Polybrominated biphenyls (PBB), Polybrominated diphenyl ethers (PBDE), Bis(2-ethylhexyl) phthalate (DEHP), Butyl benzyl phthalate (BBP), Dibutyl phthalate (DBP), Diisobutyl phthalate (DIBP).

In the case of individual assemblies, maximum concentrations maybe exceeded within the permissible exceptions of the Directive.

Schwabach, 04. September 2024

Andreas Hahn
Managing Director

Jörg Ziel
Quality Manager

China RoHS DECLARATION OF CONFORMITY

Heidolph Scientific Products GmbH has made reasonable efforts to ensure that hazardous materials and substances may not be used in its products.

In order to determine the concentration of hazardous substances in all homogeneous materials of the subassemblies, a "Product Conformity Assessment" (PCA) procedure was performed. As defined in GB/T 26572 the "Maximum Concentration Value" limits (MCV) apply to these restricted substances:

- Lead (Pb): 0.1%
- Mercury (Hg): 0.1%
- Cadmium (Cd): 0.01%
- Hexavalent chromium (Cr(VI)): 0.1%
- Polybrominated biphenyls (PBB): 0.1%
- Polybrominated diphenyl ether (PBDE): 0.1%

Environmental Friendly Use Period (EFUP)

EFUP defines the period in years during which the hazardous substances contained in electrical and electronic products will not leak or mutate under normal operating conditions. During normal use by the user such electrical and electronic products will not result in serious environmental pollution, cause serious bodily injury or damage to the user's assets.

The Environmental Friendly Use Period for Heidolph Scientific Products GmbH products is 25 years.

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This table is created according to SJ/T 11364-2014

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Note: Table applies to all products. Some of the components or parts listed above may not be part of the enclosed product.

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O: Indicates that the above mentioned hazardous substance contained in all homogeneous materials of the part is below the required limit as defined in GB/T 26572.

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X: Indicates that the above mentioned hazardous substance contained in at least one of the homogeneous materials of this part is above the required limit as defined in GB/T 26572.

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Apart from the disclosures in the above table, the subassemblies are not intentionally manufactured or formulated with lead (Pb), mercury (Hg), cadmium (Cd), hexavalent chromium (CrVI), polybrominated biphenyls (PBB), and polybrominated diphenyl ethers (PBDE).

Products manufactured by Heidolph Scientific Products GmbH may enter into further devices or can be used together with other appliances.

With these products and appliances in particular, Heidolph Scientific Products GmbH will not take responsibility for the EFUP of those products and appliances.

Schwabach, 04. September. 2024

Andreas Hahn
Managing Director

Jörg Ziel
Quality Manager

© Heidolph Scientific Products GmbH

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Subject to change without notice. The printed version of this document is not regularly updated. The latest issue of this document can be found by visiting our homepage.

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MATERIAL CONTENT DECLARATION FOR Heidolph Scientific Products GmbH PRODUCTS