# Temperature

Controlled Shaking

Betriebsanleitung  
Operating instructions  
Notice d'instructions

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\_\_\_\_\_ Datum

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\_\_\_\_\_ Firmenstempel

Translation of the original instructions

Page 24 – 39

Traduction de la notice originale

Page 44 – 59

Zertifikate/Certifications

## Contents

About this document .....	24
Typographic conventions .....	24
Copyright protection .....	24
Safety instructions in the official languages of the European Union .....	24
Directives applied, product certification .....	25
Copyright notice .....	25
California Residents .....	25
Residual risk .....	25
Intended use .....	25
Compliant use .....	25
Reasonably foreseeable misuse .....	26
Transportation .....	26
Storage .....	26
Acclimatization .....	26
Permissible ambient conditions .....	26
Energy saving potential and cost efficiency .....	26
General safety instructions .....	27
Electrical safety .....	27
Data security .....	27
Operational safety .....	27
Occupational safety .....	28
Personal protective equipment (PPE) .....	28
Environmental protection .....	28
Biohazard .....	28
Other regulations .....	28
Mechanical design .....	29
Positioning the device .....	30
Power supply .....	30
Mount/dismantle test tube tray and accessories .....	31
Switching the device on/off .....	31
Shaking modes .....	32
Automatic mode .....	32
Continuous operation .....	32
Regulate shaking movement intensity .....	33

Troubleshooting.....	34
Technical specifications .....	35
Scope of delivery .....	36
Accessories .....	36
Device service.....	37
General cleaning instructions.....	37
Repairs - Return of equipment.....	37
Maintenance.....	38
Disposal.....	38
Warranty statement.....	38
Contact details .....	38
Certificate of decontamination.....	39

## About this document

This operating manual describes the features and operation of test tube shakers type Hei-MIX Reax Top and Hei-MIX Reax Control. The operating instructions are an integral part of the described device!

## Typographic conventions

This document uses the following symbols, signal words, and highlights:

	Description
	<p>Warning symbols in combination with a signal word indicate dangers:</p> <p><b>DANGER</b> Indicates an immediate dangerous situation. Failure to respect the indications will result in death or serious injury.</p> <p><b>WARNING</b> Indicates a potential danger. Failure to respect the indications will result in serious injuries.</p> <p><b>CAUTION</b> Indicates a potential hazard. If not avoided, damage to property and minor to moderate injuries can occur.</p>
	<p>Mandatory signs are used to indicate important and useful information on handling a product. This information is used to ensure operational safety and to maintain the value of the product.</p>
[ ... ]	<p>Square brackets identify labels of control elements on the device as well as labels and entries in software masks and user interfaces.</p>
→	<p>The arrow indicates specific instructions to be followed to ensure operational safety when handling the product.</p>

## Copyright protection

This document is protected by copyright and is intended for use by the purchaser of the product only.

No transfer to third parties, reproduction in any form, including excerpts, and by any means, as well as utilization and/or disclosure of the contents is permitted without the prior written consent of Heidolph Scientific Products GmbH. Any violation is subject to compensation for damage.

## Safety instructions in the official languages of the European Union

A summary of all safety instructions in the official languages of the European Union can be found in our Safety Guide for the product group Hei-SHAKE (shakers and mixers, Ref. 01-005-006-99). This document is available for download on our homepage in the most up-to-date version.

## Directives applied, product certification

### CE marking

The device meets all requirements of the following directives:

- Machinery Directive 2006/42/EG
- EMC Directive 2014/30/EU

## Copyright notice

The software implemented in this product is protected by copyright laws. The rights holder is Heidolph Scientific Products GmbH, Walpersdorfer Straße 12, 91126 Schwabach/Germany. Any open source software components in this software are excluded from our copyright. Further information is available in the service area on our website [www.heidolph.com](http://www.heidolph.com).

Violations of copyright (e.g. unauthorized use or modification of the software) may result in civil claims (e.g. omission, compensation for damages) and/or entail criminal penalties within the meaning of the legislation of the destination country.

## California Residents

Important information for California residents regarding Prop 65:  
Please visit [www.P65Warnings.ca.gov](http://www.P65Warnings.ca.gov) for more information.

## Residual risk

The device has been designed and manufactured in accordance with the state-of-the-art standards at the time of development and the recognized safety regulations. During mounting and use, as well as during maintenance, repair and cleaning work, there are nevertheless certain residual risks associated with the described device.

These are identified and described at the appropriate points in this document.

## Intended use

The device has been developed and designed by the manufacturer for the following laboratory activities:

- Shaking
- Mixing
- Emulsifying
- Suspending
- Dissolving

The areas of application of the described platform shaker include chemical, biological and environmental-analytical laboratory and research applications, applications of basic research and comparable facilities.

Due to its design, the device in its delivery condition may only be used in analytical processes or in laboratory-like conditions in the food, cosmetics, and pharmaceutical industries as well as other comparable industries that manufacture products intended for consumption by humans or animals, or for use on humans or animals.

Any other use of this device is not considered as intended!

## Compliant use

The compliance of each individual application must always be evaluated by the user. Any additional measures necessary to ensure compliance are always the responsibility of the user.

## Reasonably foreseeable misuse

Additional measures may be necessary for use under conditions or for purposes deviating from the intended use, and/or specific guidelines and safety regulations must be observed. Corresponding requirements must be evaluated and implemented by the operator in each individual case.

Compliance with and implementation of all relevant directives and safety measures for the respective field of application is the sole responsibility of the operator.

All risks resulting from improper use are borne solely by the operator.

The device may only be operated by authorized and instructed personnel. Training and qualification of the operating personnel as well as ensuring that the device is handled responsibly are the sole responsibility of the operator!

## Transportation

During transportation, avoid severe shocks and mechanical stresses that can cause damage to the device. Keep the original packaging in a dry and protected place for later use.

## Storage

Always store the product in its original packaging. To protect against damage and excessive material aging, store the device in an environment that is as dry, temperature-stable and dust-free as possible.

## Acclimatization

After each transportation and after storage under critical climatic conditions (e.g., high temperature difference between inside and outside), allow the device to acclimatize at room temperature for at least two hours to prevent possible damage due to condensation before commissioning it in the place of use. If necessary, extend the acclimatization phase if the temperature differences are very high.

Make all supply connections only after the device has been acclimatized!

## Permissible ambient conditions

The device is designed for indoor use only. The device is **NOT** suitable for outdoor use! The device is **NOT** suitable for use in potentially explosive areas!

When used in corrosive atmospheres, the service life of the device may be reduced depending on the concentration, duration and frequency of exposure.

## Energy saving potential and cost efficiency

Even in standby mode, small amounts of energy are used to supply individual device functions. However, especially over longer periods of nonuse, this resource consumption adds up and causes avoidable costs.

Therefore, always switch off the device when not in use, especially for long periods of time (devices with main switch: Main switch in position [O]; devices without main switch: Unplug) to effectively avoid unnecessary energy consumption and the resulting costs.

## General safety instructions

Before commissioning and using the device, familiarize yourself with all the safety regulations and occupational safety guidelines applicable at the place of use and observe them at all times.

Only operate the device if it is in faultless technical condition. In particular, ensure that there is no visible damage on the device itself and, where applicable, on connected devices or the supply connections.

If there is missing or misleading information on the device or regarding occupational safety, contact the responsible safety specialist or our technical service.

Only use the device in accordance with the regulations for intended use.

## Electrical safety

Before connecting the device to the power supply, ensure that the voltage indicated on the rating plate matches the specifications of the local power utility company.

Ensure that the power supply circuit provided is protected by means of a residual-current device (RCD).

Always use the power supply cord provided with the device.

Prior to use, check that the device and the power supply cord are free of visible damage.

Make sure that the power plug can be reached directly at any time, in order to disconnect the device from the power supply at any time without delay.

Have repairs and/or maintenance work on the device carried out exclusively by an authorized electrician or by the technical service department of Heidolph Scientific Products GmbH.

Always disconnect the device from the power supply system before carrying out any maintenance, cleaning or repair work.

## Data security

The user is responsible for ensuring data security when transferring data between the described device and other devices.

Only use secure networks for the data transfer and avoid use of critical infrastructure.

Only use high-quality shielded data cables for the data transfer.

For data transfer via USB, an industrial standard USB hub should be preferably used to ensure the most stable connection possible.

## Operational safety

Operate the device under a closed ventilated fume hood when working with potentially hazardous substances (see EN 14175 and DIN 12924).

Do not make any unauthorized changes or modifications to the device!

Only use genuine spare parts and accessories, or those expressly approved by the manufacturer!

Rectify malfunctions or faults on the device immediately.

Switch off and disconnect the device from the power supply, preventing reconnection, if it is not possible to eliminate the malfunction or rectify the fault immediately.

Observe all other applicable regulations such as laboratory and workplace guidelines, recognized safety technology rules and special local regulations.

## Occupational safety

Always use the prescribed personal protective equipment (PPE) such as protective clothing, safety goggles, protective gloves, safety shoes, etc.

Do not operate any other devices in the immediate vicinity of the device ...

- which can generate electromagnetic fields in the frequency range between  $9 \times 10^3$  Hz to  $3 \times 10^{11}$  Hz,
- which generate emission or radiation sources in the frequency range  $3 \times 10^{11}$  Hz to  $3 \times 10^{15}$  Hz (in the optical spectral range wavelengths from 1,000  $\mu\text{m}$  to 0.1  $\mu\text{m}$ ),
- which generate ultrasonic or ionizing waves.

Do not process any substances that could release energy in an uncontrolled manner (e.g. self-ignition).

Do not process substances in which the energy input through mixing poses a danger.

Wipe off any liquid that may have spilled on the device immediately.

## Personal protective equipment (PPE)

The operating company must determine and provide the necessary PPE, depending on the respective application and the media and chemicals used.

The required instruction of the personnel is solely within the operating company's responsibility.

## Environmental protection

When processing environmentally hazardous substances, take appropriate measures to avoid hazards to the environment.

The evaluation of appropriate measures such as the marking of a hazardous area, their implementation, and the training of the relevant personnel is the sole responsibility of the operator!

## Biohazard

When processing biohazardous substances, take appropriate measures to prevent hazards to persons and the environment, including:

- Instruction of the personnel regarding the necessary safety measures.
- Provision of personal protective equipment (PPE) and instruction of the personnel in its use.
- Marking of the device with the biohazard warning symbol.

The evaluation of appropriate measures such as the marking of a hazardous area, their implementation, and the training of the relevant personnel is the sole responsibility of the operator!

## Other regulations

In addition to the notes and instructions in this document, observe all other applicable regulations such as laboratory and workplace guidelines, hazardous substances ordinances, recognized rules of safety engineering and occupational medicine as well as particular local regulations!

Noncompliance will invalidate any warranty claims against Heidolph Scientific Products GmbH.

The operating company is solely liable for all damage resulting from unauthorized changes or modifications to the device, from the use of unapproved or non-genuine spare parts and accessories, or from disregarding the safety instructions and hazard warnings or the manufacturer's instructions!

## Mechanical design

### Hei-MIX Reax Top

- 1 Mounting thread for tube holder
- 2 Test tube tray (max.  $\varnothing$ : 20 mm)
- Slide switch: change between automatic and continuous operation
- 4 On/off switch
- 5 Speed controller

### Hei-MIX Reax Control

- 1 Mounting thread for tube holder
- 2 Test tube tray (max.  $\varnothing$ : 20 mm)
- Slide switch: change between automatic and continuous operation
- 4 On/off switch
- 5 Speed controller

### Positioning the device

---

#### **CAUTION**

Improper positioning of the device may result in direct and indirect damage to property due to falling and/or spilling of fluids!

Observe the instructions for the correct positioning of the device.

Observe the generally valid rules for the safe positioning of work equipment.

---

- The surface must be even, non-slip, proper and sufficiently stable.
- The device must be freely accessible at all times.
- Make sure that a sufficient safety distance to moving components can be maintained on all sides.

The professional mounting and positioning of the device including the supplied accessories is within the sole responsibility of the operator!

Heidolph Scientific Products GmbH shall under no circumstances be liable for any personal injury or damage to property resulting from failure to observe the instructions for the correct installation and positioning of the device.

### Power supply

---

#### **DANGER**

Observe the notes and instructions in the section on "Elektrische Sicherheit" on page 7.

---

#### **Connect the power supply cord**

- First, connect the supplied power supply cord to the appliance plug.
- Then connect the power plug to a grounded mains socket-outlet.

#### **Disconnect the mains connection cord**

- First, remove the mains plug from the mains socket-outlet.
- Then disconnect the power supply cord from the device connector.

### Mount/dismantle test tube tray and accessories

The test tube tray as well as the available accessories (see “Accessories” on page 36) can be removed and attached without tools of any kind. The following description refers by way of example to the test tube tray supplied as standard. Proceed in the same way for the accessories.

#### Removing the test tube tray

- Grasp the soft edge of the test tube tray and pull it up.

#### Mounting the test tube tray

- When placing the test tube tray, make sure that the groove slides smoothly over the tongue on the device.
- Push the test tube tray onto the device with light pressure until it stops.

#### Switching the device on/off

- Set the speed to zero before switching on the device. To do this, turn the speed controller fully anti-clockwise.
- Use the on/off switch to turn the device on or off.
  - When on, the on/off switch lights up green.

## Shaking modes

The device can be used in automatic or continuous mode:

- In automatic mode, shaking begins as soon as a suitable test tube is pressed into the test tube tray.
- In continuous mode, the movement starts without applying pressure on the test tube tray. The shaking movement must be stopped manually!

### WARNING

In continuous mode, the device automatically restarts after an interruption (e.g. fuse tripping, power failure). There is a risk of injury in the area of movement of the device!

In the event of a power failure, switch off the device and keep sufficient distance!

### CAUTION

Depending on the load and configuration, depending on the speed, there is a risk of excessive vibration of the device during operation. There is a risk that the device moves uncontrollably, falls or tips over.

Observe the instructions for the correct positioning of the device!

Set the speed controller to zero before switching on the device. Gradually increase the speed to the desired value as the shaking movement starts. Observe the device for sufficient stability.

Check and correct the configuration if excessive vibrations are detected!

## Automatic mode

Switching between automatic and continuous mode is done via the slide switch of the device, see "Mechanical design" on page 29.

Move the slide switch [1] to the right into the [automatic mode] position, see figure. The automatic mode icon [2] must be visible:

- Place the test tube in the test tube tray and hold it firmly with a slight downward pressure.
  - The shaking movement starts by applying pressure with the test tube.
- As soon as the test tube is removed from the test tube tray, the shaking movement will stop.

## Continuous operation

Move the slide switch [1] to the left into the [continuous mode] position, see figure. The continuous mode icon [2] must be visible:

- The shaking movement starts as soon as the slide switch is in the [continuous mode] position!
- Move the slide switch to the right into the [automatic mode] position or switch off the device to stop the movement.

## Regulate shaking movement intensity

The shaking movement intensity is set by the speed and by means of the speed controller of the device.

### WARNING

If the shaking operation starts too quickly, fluid may splash out of the test tube.

Use recommended personal protective equipment (e.g. safety glasses, protective clothing).

Set the speed controller to zero before switching on the device. Gradually increase the speed to the desired value as the shaking movement starts.

### WARNING

Depending on the effort and duration, holding the sample in automatic mode can lead to overloads in the hand and arm area.

When using the device, observe the specifications of the EC directive 2002-44 on the permissible vibration load.

### Reax Top

- Turn the speed controller clockwise to increase the shaking movement intensity.
- Turn the speed controller counter-clockwise to decrease the shaking movement intensity.

The Reax Top model is equipped with an analogue scale with reference points:

### Reax Control

- Turn the speed controller clockwise to increase the shaking movement intensity.
- Turn the speed controller counter-clockwise to decrease the shaking movement intensity.

The Reax Control model is equipped with a graduated scale:

## Troubleshooting

Malfunction	Possible cause	Remedy
On/off switch remains dark after switching on	No power supply	<ul style="list-style-type: none"> <li>→ Check that the power plug is correctly connected</li> <li>→ Check circuit breaker</li> </ul>
	LED defective	<ul style="list-style-type: none"> <li>→ Contact authorized dealer or technical service</li> </ul>
Shaking movement does not start in automatic mode	Slide switch not in correct position	<ul style="list-style-type: none"> <li>→ Push the slide switch all the way to the right</li> </ul>
	Not enough pressure on the test tube tray	<ul style="list-style-type: none"> <li>→ Apply more pressure with test tube</li> </ul>
	Too much pressure on the test tube tray, device is switched off within 60 s	<ul style="list-style-type: none"> <li>→ Reduce pressure on the test tube tray</li> </ul>
	Test tube too slanted	<ul style="list-style-type: none"> <li>→ Center test tube on tray</li> <li>→ If possible, switch to continuous mode</li> </ul>
Acoustic signal (Reax Control only)	Power recovery after interruption in continuous mode	<ul style="list-style-type: none"> <li>→ The signal tone is switched off after approx. 2 seconds, shaking movement is continued</li> </ul>
Shaking movement stops in continuous mode	<ul style="list-style-type: none"> <li>→ Motor defective</li> <li>→ Switch defective</li> <li>→ Electronics defective</li> </ul>	<ul style="list-style-type: none"> <li>→ Contact authorized dealer or technical service</li> </ul>
	<ul style="list-style-type: none"> <li>→ Overheat protection triggered</li> </ul>	<ul style="list-style-type: none"> <li>→ Allow the device to cool, restart, reduce load</li> </ul>

If a fault cannot be rectified with the described suggestions, please contact an authorized sales representative or our technical service (see section "Contact details" on page 38).

## Technical specifications

Variants	Reax Top 230 V 50 Hz	100 – 2,500 rpm
	Reax Top 230 V 60 Hz	200 – 3,200 rpm
	Reax Top 115 V 60 Hz	
	Reax Control 230 V 50/60 Hz	0 – 2,500 rpm
Reax Control 115 V 50/60 Hz		
Power input	51 W	
Protection class (IEC 61140)	1 Ⓢ	
Protection class (IEC 60529)	IP 22	
Acoustic pressure (dB(A)) (in accordance with IEC 61010)	< 50 to 2,500 rpm < 58 to 3,200 rpm	
Drive	Shaded pole motor	
Overheat protection	self-resetting	
Motion	circular vibrating	
Orbit	5 mm	
Operating mode	Automatic and continuous operation	
Weight	2.8 kg	
Dimensions (W x H x D)	134 x 105 x 172 mm	
<b>Hei-Mix model</b>	<b>Reax Top</b>	<b>Reax Control</b>
Speed control	analogue, scale of reference points	analogue, graduated scale
Control accuracy	unregulated	±2 %

## Scope of delivery

Component / Variants		Quantity	Product number
Hei-MIX	Reax Top*	1	541-10000-00
	Reax Control*	1	541-11000-00
Test tube tray, small		1	23-07-06-05-01
Power supply cord		1	country specific

\* the product numbers shown are for 230 V EU devices. For product numbers of variants, please contact Heidolph Scientific Products GmbH.

## Accessories

Component / Variant	Quantity	Item number
Test tube tray, large	1	549-19000-00
Support rod with vessel holder	1	549-20000-00
Holding rack for 10 test tubes	1	549-01000-00
Test tube stand (max. 6 Eppendorf vessels)	1	594-04000-00

Further accessories can be found on our homepage at [www.heidolph.com](http://www.heidolph.com)

## Device service

When carrying out service work on the device (cleaning, maintenance, repair), observe the general instructions and safety information described in this section.

### DANGER

Switch the device's main switch off and disconnect it from the power supply before carrying out maintenance work, cleaning, or repairs.

When cleaning, avoid the penetration of liquids.

Before replacing the fuses, switch off the device and disconnect the power supply cord.

Always replace the two device fuses in pairs with original manufacturer fuses. Further information on available accessories can be found on our website at [www.heidolph.com](http://www.heidolph.com)!

After fuse replacement, check the device for a safe condition according to IEC 61010-1.

## General cleaning instructions

Wipe all surfaces and the control panel with a damp cloth if necessary. Persistent contamination can be removed with mild soapy water.

### CAUTION

Clean the device's surfaces with a soft, lint-free and only slightly moistened cloth.

Never use any aggressive or abrasive cleaning agents or aids.

## Repairs - Return of equipment

Repairs to the device may only be carried out by authorized experts! Unauthorized repairs during the warranty period will result in the loss of the warranty claim. Regardless of the warranty claim, the owner is solely liable for damage caused by unauthorized repairs.

- In case of repair and before returning your device, contact our technical service at the following e-mail address:
  - [service@heidolph.de](mailto:service@heidolph.de).
- In your message, please provide us with the following information in addition to an error description:
  - Item number
  - Serial no.

The required data can be found on the rating plate of the device.

A service representative will contact you as soon as possible to agree on the next steps.

Include the completed certificate of decontamination with every device return, see "Unbedenklichkeitserklärung" on page 28.

## Maintenance

The device contains no user-serviceable components. If necessary (in the event of abnormal operating behavior such as excessive noise or heat generation, for example) contact our technical service.

## Disposal

- When disposing of the device, observe the provisions of the WEEE Directive 2012/19/EU and its transposition into national law in the country of use.
- When disposing of portable batteries, observe the provisions of the European Battery Directive 2013/56/EU and their implementation in national law in the country of use.
- Check the device and all components for residues of substances that are hazardous to health, the environment and biohazardous before disposal.
- Properly remove and dispose of residues of substances that are hazardous to health, the environment and biohazardous!

### Registration in accordance with the Electrical and Electronic Equipment Act (ElektroG)

Heidolph Scientific Products GmbH, headquartered in 91126 Schwabach/Germany, address Walpersdorfer Str. 12, is registered under the number DE 50705753 in the electronic waste equipment register of the foundation ear ([www.ear-system.de](http://www.ear-system.de)).

## Warranty statement

Heidolph Scientific Products GmbH provides a three-year warranty against material and manufacturing defects.

Glass and wear parts, transportation damage, and damage resulting from improper handling or non-intended use of the product are excluded from the warranty.

The warranty period for registered products begins on the date of purchase. Register the product with the enclosed warranty card or on our homepage [www.heidolph.com](http://www.heidolph.com).

For non-registered products, the warranty period begins with the date of the serial production (to be determined by the serial number).

In the event of material or manufacturing defects, the product will either be repaired or replaced free of charge within the warranty period.

## Contact details

### Heidolph Scientific Products GmbH

Technical service  
Walpersdorfer Str. 12  
D-91126 Schwabach/Germany  
Email: [service@heidolph.de](mailto:service@heidolph.de)

### Representations

To find your local Heidolph distributor please visit [www.heidolph.com](http://www.heidolph.com)

## Certificate of decontamination

Enclose the certificate of decontamination, duly completed, with your device return. Submissions without a certificate of decontamination cannot be processed!

# CERTIFICATE OF DECONTAMINATION

IN CASE OF RETURNS

Please fill in the required fields.

**Note: The sender must package the goods properly and appropriately for transport.**

Heidolph Scientific Products GmbH

Walpersdorfer Straße 12  
91126 Schwabach

Phone: +49 (0) 9122 9920-380

Fax: +49 (0) 9122 9920-19

E-Mail: service@heidolph.de

### SENDER

Name \_\_\_\_\_ First name \_\_\_\_\_  
 Company/institution \_\_\_\_\_ Department \_\_\_\_\_  
 \_\_\_\_\_ Workgroup \_\_\_\_\_  
 Address \_\_\_\_\_  
 ZC/City \_\_\_\_\_  
 Country \_\_\_\_\_ Phone \_\_\_\_\_  
 Email \_\_\_\_\_

### DEVICE DETAILS

Article number \_\_\_\_\_ Serial no. \_\_\_\_\_

Ticket number \_\_\_\_\_

Reason for sending in \_\_\_\_\_  
 \_\_\_\_\_

Has the device been cleaned, decontaminated/disinfected? **Yes** **No** (Please mark as applicable)

If yes, which measures were carried out?  
 \_\_\_\_\_

Does this device pose a risk to people and/or the environment due to the processing of substances that are hazardous to health, the environment and/or are biohazardous? **Yes** **No** (Please mark as applicable)

If yes, with which substances did the device come into contact?  
 \_\_\_\_\_

### LEGALLY BINDING DECLARATION

The principal/consignor is aware that they are liable to the agent/consignee for losses or damage incurred due to incomplete and incorrect information.

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hereby declare, that the product (from serial number on 200066855) designated below is in compliance with the basic requirements of all applicable EU-directives stated below with regard to design, type of model sold and manufactured by us. This certificate will be invalid if the product is modified without the prior written consent and agreement of the manufacturer.

Reax 2	541-21001-XX		
Reax top	541-10000-XX		
Reax control	541-11000-XX		
Reax 20/4	541-20004-XX		
Reax 20/8	541-20008-XX		
Reax 20/12	541-20012-XX		
Titramax 100	544-11100-XX, 544-11200-XX		
Duomax 1030	543-32105-XX, 543-32205-XX, 543-32210-XX		
Polymax 1040	543-42105-XX, 543-42110-XX, 543-42205-XX, 543-42210-XX		
Polymax 2040	542-40005-XX, 542-40010-XX		
Multi Reax	545-10000-XX		
Vibramax 100	544-21200-XX	Titramax 101	544-11300-XX
Rotamax 120	544-41200-XX	Vibramax 110	544-31200-XX
Titramax 1000	544-12200-XX	Inkubator 1000	549-90010-XX
Unimax 1010	543-12310-XX	Promax 1020	543-22332-XX
Unimax 2010	542-10020-XX	Promax 2020	542-20020-XX
Hei SHAKE Orbital Core	546-11XXX-XX		

Maschinenrichtlinie / Machinery Directive 2006/42/EG  
EMV-Richtlinie / Electromagnetic Compatibility Directive 2014/30/EU  
Delegierte (RoHS-) Richtlinie / Delegated (RoHS) Directive 2015/863/EU  
Angewandte (harmonisierte) Normen / (Harmonized) Standards applied:  
EN ISO 12100:2010, EN 61326-1:2013, EN 61010-1:2010, EN IEC 63000:2018

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**UK  
CA**

## Declaration of Conformity

In accordance with UK Government guidance

This declaration is issued under the sole responsibility of the manufacturer,

**Heidolph Scientific Products GmbH**

Walpersdorfer Straße 12

91126 Schwabach / Germany

Product: Shaker & Mixer

Model:

Reax top /control	541-1XXXX-XX
Reax 20/4...12	541-200XX-XX
Reax 2	541-21001-XX
Duomax 1030	543-32XXX-XX
Unimax X010	54X-1XXX0-XX
Promax X020	54X-2XXXX-XX
Polymax X040	54X-4XXXX-XX
Titramax 1XXX	544-1XXXX-XX
Vibramax 1XX	544-X1200-XX
Rotamax 120	544-41200-XX
Multi Reax	545-10000-XX
Inkubator 1000	549-90010-XX
Hei SHAKE Orbital Core	546-11XXX-XX

Description:

Shaker & Mixer with different working directions

The object of the declaration described above is in conformity with the relevant UK Statutory Instruments (and their amendments):

2008 No. 1597

2008 2016 No. 1091

2012 No. 3032

*The Supply of Machinery (Safety) Regulations*

*The Electromagnetic Compatibility Regulations 2016*

*The Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic Equipment Regulations 2012*

and complies with the following technical standards :

EN ISO 12100:2010, EN 61326-1:2013, EN 61010-1:2010, EN IEC 63000:2018

UK Authorised Representative (for authorities only)

ProductIP ( UK ) Ltd.

8. Northumberland Av.

London WC2N 5BY

Signed for and on behalf of Heidolph Scientific Products GmbH

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With this declaration, we confirm (according to current knowledge) that all sold laboratory devices and components by Heidolph Scientific Products GmbH fulfill the requirements of the EU directive 2011/65/EU (RoHS) and the extension 2015/863. All devices are compatible with the requirements of the RoHS Directive for the following materials:

Max. 0,01% of weight in homogenous material for cadmium and max. 0,1% of the weight in homogeneous material for Lead, Mercury, Hexavalent chromium, Polybrominated biphenyls (PBB), Polybrominated diphenyl ethers (PBDE), Bis(2-ethylhexyl) phthalate (DEHP), Butyl benzyl phthalate (BBP), Dibutyl phthalate (DBP), Diisobutyl phthalate (DIBP).

In the case of individual assemblies, maximum concentrations maybe exceeded within the permissible exceptions of the Directive.

Schwabach, 04. September 2024

Handwritten signature of Andreas Hahn in blue ink.

Andreas Hahn  
Managing Director

Handwritten signature of Jörg Ziel in blue ink.

Jörg Ziel  
Quality Manager

## China RoHS DECLARATION OF CONFORMITY

Heidolph Scientific Products GmbH has made reasonable efforts to ensure that hazardous materials and substances may not be used in its products.

In order to determine the concentration of hazardous substances in all homogeneous materials of the subassemblies, a "Product Conformity Assessment" (PCA) procedure was performed. As defined in GB/T 26572 the "Maximum Concentration Value" limits (MCV) apply to these restricted substances:

- Lead (Pb): 0.1%
- Mercury (Hg): 0.1%
- Cadmium (Cd): 0.01%
- Hexavalent chromium (Cr(VI)): 0.1%
- Polybrominated biphenyls (PBB): 0.1%
- Polybrominated diphenyl ether (PBDE): 0.1%

### Environmental Friendly Use Period (EFUP)

EFUP defines the period in years during which the hazardous substances contained in electrical and electronic products will not leak or mutate under normal operating conditions. During normal use by the user such electrical and electronic products will not result in serious environmental pollution, cause serious bodily injury or damage to the user's assets.

The Environmental Friendly Use Period for Heidolph Scientific Products GmbH products is 25 years.

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This table is created according to SJ/T 11364-2014

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**Note:** Table applies to all products. Some of the components or parts listed above may not be part of the enclosed product.

- O: 表示该有毒有害物质在该部件所有均质材料中的含量均在GB/T 26572规定的限量要求以下。
- O: Indicates that the above mentioned hazardous substance contained in all homogeneous materials of the part is below the required limit as defined in GB/T 26572.
- X: 表示该有毒有害物质至少在该部件某一均质材料中的含量超出GB/T 26572规定的限量要求。
- X: Indicates that the above mentioned hazardous substance contained in at least one of the homogeneous materials of this part is above the required limit as defined in GB/T 26572.

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Apart from the disclosures in the above table, the subassemblies are not intentionally manufactured or formulated with lead (Pb), mercury (Hg), cadmium (Cd), hexavalent chromium (CrVI), polybrominated biphenyls (PBB), and polybrominated diphenyl ethers (PBDE).

Products manufactured by Heidolph Scientific Products GmbH may enter into further devices or can be used together with other appliances .

With these products and appliances in particular, Heidolph Scientific Products GmbH will not take responsibility for the EFUP of those products and appliances.

Schwabach, 04. September. 2024

Andreas Hahn  
Managing Director

Jörg Ziel  
Quality Manager

© Heidolph Scientific Products GmbH

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Subject to change without notice. The printed version of this document is not regularly updated. The latest issue of this document can be found by visiting our homepage.

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MATERIAL CONTENT DECLARATION FOR Heidolph Scientific Products GmbH PRODUCTS