
# Precise

Dosing and Suspending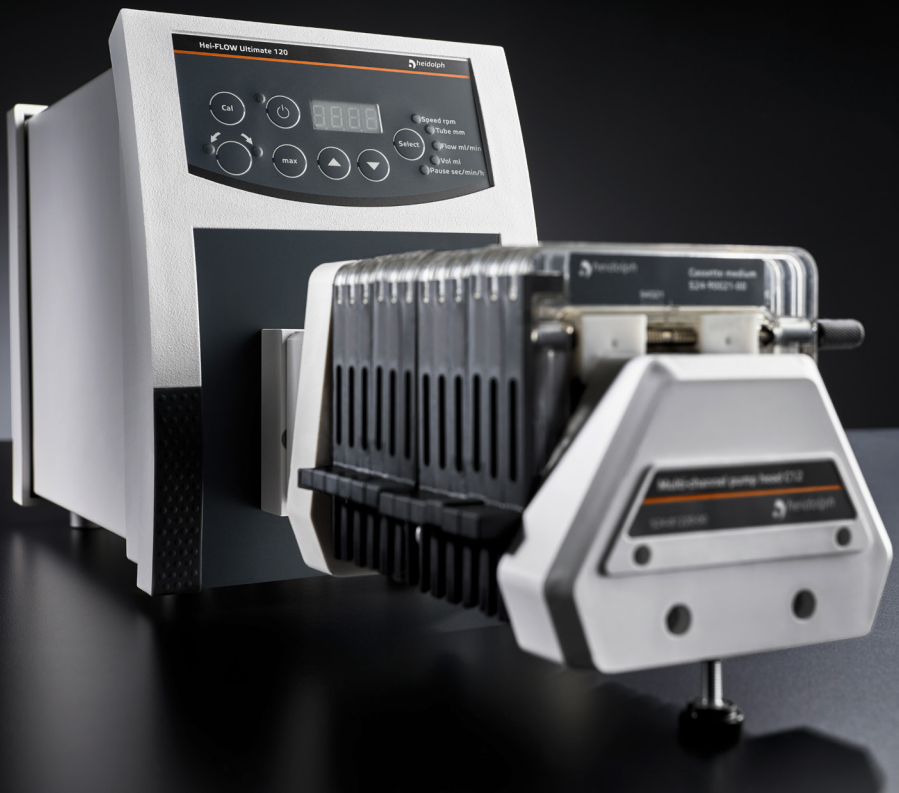

## Safety Guide

Peristaltikpumpen  
Peristaltic pumps  
Pompes péristaltiques

## Index

German .....	3	
English .....	8	
Български .....	12	
Čeština .....	17	
Dansk .....	21	
Eesti .....	25	
Español .....	29	
Suomi .....	33	
French .....	37	
Ελληνικά .....	42	
Hrvatski .....	47	
Magyar .....	51	
Italiano .....	55	
Lietuvių kalba .....	60	
Latviešu .....	64	
Nederlands .....	68	
Polski .....	73	
Português .....	78	
Română .....	83	
Svenska .....	88	
Slovenščina .....	92	
Slovensky .....	96	

## Zu diesem Dokument

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## About this document

This document contains a summary of all safety instructions for Heidolph Scientific Products GmbH peristaltic pumps in the official languages of the European Union.

This document in no way replaces the operating instructions supplied with the product!

For more detailed information, please refer to the original instructions (in German) or the translation of the original instructions supplied with the product.

## Typographic conventions

Standardized symbols, signal words and highlighting elements are used in this document to warn of hazards and to identify important information or rather special text contents.

Symbol	Signal word / explanation
	Warning symbols in combination with a signal word indicate dangers: <b>DANGER</b> Indicates an immediate dangerous situation. Failure to respect the indications will result in death or serious injury.
	<b>WARNING</b> Indicates a potential danger. Failure to respect the indications will result in serious injuries.
	<b>CAUTION</b> Indicates a potential hazard which, if not avoided, damage to property and minor to moderate injuries can occur.
	Mandatory signs are used to indicate important and useful information on handling a product. This information is used to ensure operational safety and to maintain the value of the product.

## Copyright protection

This document is protected by copyright and is intended for use by the purchaser of the product only.

No transfer to third parties, reproduction in any form, including excerpts, and by any means, as well as utilization and/or disclosure of the contents is permitted without the prior written consent of Heidolph Scientific Products GmbH. Any violation is subject to compensation for damage.

## Intended use

Peristaltic pumps of the company Heidolph Scientific Products GmbH were especially developed for conveying and dosing liquids.

Detailed information can be seen in the associated operating instructions. The areas of application of Heidolph Scientific Products GmbH' peristaltic pumps include especially chemical, biological and environmental-analytical laboratory and research applications, applications of basic research and comparable facilities.

Due to their design, the peristaltic pumps of Heidolph Scientific Products GmbH in their delivery condition may only be used in analytical processes or in laboratory-like conditions in the food, cosmetics, and pharmaceutical industries as well as other comparable industries that manufacture products intended for consumption by humans or animals, or for use on humans or animals.

Any other use is deemed to be not as intended!

## General safety instructions

- Before commissioning and using the device, familiarize yourself with all the safety regulations and occupational safety guidelines applicable at the place of use and observe them at all times.
- Only operate the device if it is in faultless technical condition. In particular, ensure that there is no visible damage on the device itself, the power supply cord and, where applicable, on connected devices or the supply connections.
- If there is missing or misleading information on the device or regarding occupational safety, contact the responsible safety specialist or our technical service.
- Only use the device in accordance with the regulations for intended use.

### Electrical safety

- Before connecting the device to the power supply, ensure that the voltage indicated on the rating plate matches the specifications of the local power utility company.
- Ensure that the power supply circuit provided is protected by means of a residual-current device (RCD).
- Always use the power supply cord provided with the device.
- The device must only be supplied with power from a properly grounded mains socket-outlet.
- Have repairs and/or maintenance work on the device carried out exclusively by an authorized electrician or by the technical service department of Heidolph Scientific Products GmbH.
- Always disconnect the device from the power supply system before carrying out any maintenance, cleaning or repair work.

### Data security

- The user is responsible for ensuring data security when transferring data between a peristaltic pump of the company Heidolph Scientific Products GmbH and other devices/systems.
- Only use secure networks for the data transfer and avoid use of critical infrastructure.
- Only use high-quality shielded data cables for the data transfer.
- For data transfer via a USB B connection, an industrial standard USB hub should be preferably used to ensure the most stable connection possible.

### Operational safety

- Operate the device under a closed ventilated fume hood when working with potentially hazardous substances (see EN 14175 and DIN 12924).
- Do not make any unauthorized changes or modifications to the device!
- Only use genuine spare parts and accessories, or those expressly approved by the manufacturer!
- Rectify malfunctions or faults on the device immediately.
- Switch off and disconnect the device from the power supply, preventing reconnection, if it is not possible to eliminate the malfunction or rectify the fault immediately.
- Observe all other applicable regulations such as laboratory and workplace guidelines, recognized safety technology rules and special local regulations.

### Occupational safety

- Always use the prescribed personal protective equipment (PPE) such as protective clothing, safety goggles, protective gloves, safety shoes, etc.
- Persons with pacemakers or other susceptible implants must maintain a minimum distance of ten cm from the device to prevent possible interference.
- Do not operate any other devices in the immediate vicinity of the device
  - which generate emission or radiation sources in the frequency range  $3 \times 10^{11}$  Hz to  $3 \times 10^{15}$  Hz (in the optical spectral range wavelengths from 1,000  $\mu\text{m}$  to 0,1  $\mu\text{m}$ ),
  - which generate ultrasonic or ionizing waves.
- Do not process any substances that could release energy in an uncontrolled manner (e.g. self-ignition).
- Do not process substances in which the energy input through mixing poses a danger.
- Do not work with an open flame in the immediate vicinity of the device (risk of explosion or fire when processing explosive/flammable media).
- Wipe off any liquid that may have spilled on the device immediately.
- Always switch off the device or disconnect the device from the power supply when it is not in use.

### Personal protective equipment (PPE)

The operating company must determine and provide the necessary PPE, depending on the respective application and the media and chemicals used.

The corresponding instruction of the personnel is solely within the operating company's responsibility.

### Environmental protection

When processing environmentally hazardous substances, take appropriate measures to avoid hazards to the environment.

The evaluation of corresponding measures such as the marking of a hazardous area, their implementation, and the training of the relevant personnel is the sole responsibility of the operator!

### Biohazard

When processing biohazardous substances, take appropriate measures to prevent hazards to persons and the environment, including:

- Instruction of the personnel regarding the necessary safety measures.
- Provision of personal protective equipment (PPE) and instruction of the personnel in its use.
- Marking the device with a biohazard warning symbol.

The evaluation of corresponding measures such as the marking of a hazardous area, their implementation, and the training of the relevant personnel is the sole responsibility of the operator!

### Other regulations

In addition to the notes and instructions in this document, observe all other applicable regulations such as laboratory and workplace guidelines, hazardous substances ordinances, recognized rules of safety engineering and occupational medicine as well as particular local regulations!

Noncompliance will invalidate any warranty claims against Heidolph Scientific Products GmbH.

The operator is solely liable for all damage resulting from unauthorized changes or modifications to the device, from the use of unauthorized or non-genuine spare parts and accessories, or from disregarding the safety instructions and hazard warnings or the manufacturer's instructions!

## Device-specific safety instructions

### Mounting

#### CAUTION

Place the device for operation always on a clean, even and sufficiently stable work area.

The device must be freely accessible at all times.

Observe the generally valid rules for the safe installation of work equipment.

### Interfaces

#### DANGER

Penetrating liquids pose the risk of an electric shock.

In addition, the interfaces and other electronic components inside the housing can be damaged and cause the pump to fail!

- Use only the connectors recommended by the manufacturer.
- Make sure that the connection plugs are inserted correctly.
- Always protect unused interfaces with the supplied screw cap.

### Pump head mounting

#### CAUTION

When removing/mounting the pump head, observe all relevant safety regulations for mechanical activities to avoid injury to hands and fingers.

Always disconnect the device from the mains before removing/mounting the pump head to avoid unintended start-up!

Only dismantle protective and/or safety devices where necessary and exclusively for the purpose of removing/mounting components and re-assemble them properly after the work has been completed.

Before switching on the device, make sure that all protective and/or safety devices are properly mounted.

### Mounting tubes

#### CAUTION

When removing/mounting the tubes, observe all relevant safety regulations for mechanical activities to avoid injury to hands and fingers.

Always disconnect the device from the mains before installing the tube(s) to avoid injury and/or damage to property due to an unintended pump start!

Avoid excessive force when mounting the tube(s): All adjustment and fastening elements must be operated with manual force only.

Only dismantle protective and/or safety devices where necessary and exclusively for the purpose of removing/mounting components and re-assemble them properly after the work has been completed.

Before switching on the device, make sure that all protective and/or safety devices are properly mounted.

Check that all cassettes and tubes are securely seated before starting the pump.

### Device service

#### DANGER

Switch the device's main switch off and disconnect it from the power supply before carrying out maintenance work, cleaning, or repairs.

When cleaning, avoid the penetration of liquids.

Before replacing the fuses, switch off the device and disconnect the power supply cord.

Always replace the two device fuses in pairs with original manufacturer fuses. Further information on available accessories can be found on our website at [www.heidolph.com](http://www.heidolph.com)!

After fuse replacement, check the device for a safe condition according to IEC 61010-1.

#### CAUTION

Clean the device's surfaces with a soft, lint-free and only slightly moistened cloth.

Never use any aggressive or abrasive cleaning agents or aids.

### Repairs

Repairs to the device may only be carried out by authorized experts!

Unauthorized repairs during the warranty period will result in the loss of the warranty claim.

The owner is solely liable for damage caused by unauthorized repairs.

If repairs are necessary, contact an authorized dealer or our technical service.

### Maintenance

The device contains no user-serviceable components. If necessary (in the event of abnormal operating behavior such as excessive noise or heat generation, for example) contact our technical service.

### Other regulations

In addition to the notes and instructions in this document, observe all other applicable regulations such as laboratory and workplace guidelines, hazardous substances ordinances, recognized rules of safety engineering and occupational medicine as well as particular local regulations!

### Contact details Heidolph Scientific Products GmbH

#### Heidolph Scientific Products GmbH

Technical Service  
Walpersdorfer Str. 12  
D-91126 Schwabach/Germany  
Email: [service@heidolph.de](mailto:service@heidolph.de)

#### Representations

To find your local Heidolph distributor please visit [www.heidolph.com](http://www.heidolph.com)

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### Kontaktné údaje Heidolph Scientific Products GmbH
© Heidolph Scientific Products GmbH