Unlimited
Combinations

Safety Guide

Schüttel-Misch-Geräte
Shakers and Mixers
Agitateurs mélangeurs

 **heidolph**
research made easy

Index

German	2	
English	8	
Български	13	
Čeština	19	
Tysk	24	
ελληνικά	29	
Español	35	
Eesti	40	
Suomi	45	
Français	50	
Hrvatski	56	
Magyar	61	
Italiano	66	
Lietuvių k.....	71	
Latviešu	76	
Nederlands	81	
Polski	86	
Português	92	
Română	97	
Slovenčina	102	
slovensko	107	
Svenska	112	

Zu diesem Dokument

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www.heidolph.com

About this document

This document contains a summary of all safety instructions for Heidolph shakers and mixers in the official languages of the European Union.

This document in no way replaces the operating instructions supplied with the product! For more detailed information, please refer to the original instructions (in German) or the translation of the original instructions supplied with the product.

Typographic conventions

The following symbols, signal words and highlighting elements are used in this document:

Symbol	Signal word / explanation
	<p>Warning symbols in combination with a signal word indicate dangers:</p> <p>DANGER Indicates an immediate dangerous situation. Failure to respect the indications will result in death or serious injury.</p> <p>WARNING Indicates a potential danger. Failure to respect the indications will result in serious injuries.</p> <p>CAUTION Indicates a potential hazard which, if not avoided, damage to property and minor to moderate injuries can occur.</p>
	<p>Mandatory signs are used to indicate important and useful information on handling a product. This information is used to ensure operational safety and to maintain the value of the product.</p>

Copyright protection

This document is protected by copyright and is generally intended for use by the purchaser of the product. No transfer to third parties, reproduction in any form, including excerpts, and by any means, as well as utilization and/or disclosure of the contents is permitted without the prior written consent of Heidolph Scientific Products GmbH. Any violation is subject to compensation for damage.

Intended use

Shakers and mixers of the company Heidolph Scientific Products GmbH have been specially developed for the following tasks:

- Shaking
- Mixing
- Emulsifying
- Suspending
- Separating
- Dissolving
- Staining

Detailed information can be seen in the associated operating instructions. The areas of application of Heidolph Scientific Products GmbH shakers and mixers include especially chemical, biological and environmental-analytical laboratory and research applications, applications of basic research and comparable facilities.

Due to their design, the shakers and mixers of Heidolph Scientific Products GmbH in their delivery condition may only be used in analytical processes or in laboratory-like conditions in the food, cosmetics, and pharmaceutical industries as well as other comparable industries that manufacture products intended for consumption by humans or animals, or for use on humans or animals.

Any other use is deemed to be not as intended!

Compliant use

The user is generally responsible for evaluating the conformity of his application and, if necessary, for taking additional measures.

Reasonably foreseeable misuse

Additional measures may be necessary for use under conditions or for purposes deviating from the intended use, and/or specific guidelines and safety regulations must be observed. Corresponding requirements must be evaluated and implemented by the operator in each individual case.

Compliance with and implementation of all relevant directives and safety measures for the respective field of application is generally the responsibility of the operator.

All risks resulting from improper use are borne generally by the operator.

The device may generally be operated by authorized and instructed personnel. Training and qualification of the operating personnel as well as ensuring that the device is handled responsibly are the general responsibility of the operator!

Transportation

During transportation, avoid severe shocks and mechanical stresses that can cause damage to the device. Keep the original packaging in a dry and protected place for later use.

Storage

Always store the device in its original packaging. To protect against damage and excessive material aging, store the device in an environment that is as dry, temperature-stable and dust-free as possible.

Acclimatization

After each transportation and after storage under critical climatic conditions (e.g. high temperature difference between inside and outside), allow the device to acclimatize at room temperature for at least two hours to prevent possible damage due to condensation before commissioning it in the place of use. If necessary, extend the acclimatization phase if the temperature differences are very high.

Make all supply connections (power supply, tubing) only after the device has been acclimatized!

Permissible ambient conditions

The device is designed for indoor use only. The device is not suitable for outdoor use! The device is not suitable for use in potentially explosive areas!

When used in corrosive atmospheres, the service life of the device may be reduced depending on the concentration, duration and frequency of exposure.

General safety instructions

Before commissioning and using the device, familiarize yourself with all the safety regulations and occupational safety guidelines applicable at the place of use and observe them at all times.

Only operate the device if it is in faultless technical condition. In particular, ensure that there is no visible damage on the device itself and, where applicable, on connected devices or the supply connections.

If there is missing or misleading information on the device or regarding occupational safety, contact the responsible safety specialist or our technical service.

Only use the device in accordance with the regulations for intended use.

Electrical safety

Before connecting the device to the power supply, ensure that the voltage indicated on the rating plate matches the specifications of the local power utility company.

Ensure that the power supply circuit provided is protected by means of a residual-current device (RCD).

Always use the power supply cord supplied with the device. All installation work for the power supply of the device may only be carried out by an authorized electrician or by the technical service of Heidolph Scientific Products GmbH!

Prior to use, check that the device and the power supply cord are free of visible damage.

Always have repairs and/or maintenance work on the device carried out by an authorized electrician or by the technical service department of Heidolph Scientific Products GmbH.

Always switch off and disconnect the device from the power supply, preventing reconnection, before carrying out maintenance work, cleaning, or repairs.

Data security

The user is responsible for ensuring data security when transferring data between the described device and other devices.

Always use secure networks for the data transfer and avoid use of critical infrastructure.

Always use high-quality shielded data cables for the data transfer.

For data transfer via a USB B connection, an industrial standard USB hub should be preferably used to ensure the most stable connection possible.

Operational safety

Operate the device under a closed ventilated fume hood when working with potentially hazardous substances (see EN 14175 and DIN 12924).

Do not make any unauthorized changes or modifications to the device!

Always use genuine spare parts and accessories, or those expressly approved by the manufacturer!

Rectify malfunctions or faults on the device immediately.

Switch off and disconnect the device from the power supply, preventing reconnection, if it is not possible to eliminate the malfunction or rectify the fault immediately.

Observe all relevant general and safety instructions for the connected peripheral devices (observe the supplied documentation!).

Observe all other applicable regulations such as laboratory and workplace guidelines, recognized safety technology rules and special local regulations.

Occupational safety

Always use the prescribed personal protective equipment (PPE) such as protective clothing, safety goggles, protective gloves, safety shoes, etc.

Persons with pacemakers or other susceptible implants must maintain a minimum distance of ten cm from the device to prevent possible interference.

Do not operate any other devices in the immediate vicinity of the device

- which generate emission or radiation sources in the frequency range 3×10^{11} Hz to 3×10^{15} Hz (in the optical spectral range wavelengths from 1,000 μm to 0,1 μm),
- which generate ultrasonic or ionizing waves.

Do not process any substances that could release energy in an uncontrolled manner (e.g. self-ignition).

Do not process substances in which the energy input through mixing poses a danger.

Do not work with an open flame in the immediate vicinity of the device (risk of explosion or fire when processing explosive/flammable media).

Wipe off any liquid that may have spilled on the device immediately.

Always switch off the device or disconnect the device from the power supply when it is not in use.

Personal protective equipment (PPE)

The operating company must determine and provide the necessary PPE, depending on the respective application and the media and chemicals used.

The corresponding instruction of the personnel is generally within the operating company's responsibility.

Environmental protection

When processing environmentally hazardous substances, take appropriate measures to avoid hazards to the environment.

The evaluation of corresponding measures such as the marking of a hazardous area, their implementation, and the training of the relevant personnel is the general responsibility of the operator!

Biohazard

When processing biohazardous substances, take appropriate measures to prevent hazards to persons and the environment, including:

- Instruction of the personnel regarding the necessary safety measures.
- Provision of personal protective equipment (PPE) and instruction of the personnel in its use.
- Marking the device with a biohazard warning symbol.

The evaluation of corresponding measures such as the marking of a hazardous area, their implementation, and the training of the relevant personnel is the general responsibility of the operator!

Other regulations

In addition to the notes and instructions in this document, observe all other applicable regulations such as laboratory and workplace guidelines, hazardous substances ordinances, recognized rules of safety engineering and occupational medicine as well as particular local regulations!

Noncompliance will invalidate any warranty claims against Heidolph Scientific Products GmbH.

The operator is generally liable for all damage resulting from unauthorized changes or modifications to the device, from the use of unapproved or non-genuine spare parts and accessories, or from disregarding the safety instructions and hazard warnings or the manufacturer's instructions!

Device-specific safety instructions

Mounting

CAUTION

To use the device, place it on a clean, stable, level, horizontal, and heat-resistant surface.

Make sure that the power plug can be reached directly at any time, in order to disconnect the device from the power supply at any time without delay.

Before switching on, make sure that the device is sufficiently stable.

Before switching on, always make sure that all setups are centered as far as possible and are fixed on the base unit with sufficient stability.

Keep all support and contact surfaces clean and dry.

Power supply

DANGER

The device may only be supplied with the required mains voltage via a grounded mains socket-outlet (voltage see rating plate).

Only use the supplied power supply cord.

Automatic restart

WARNING

Before activating the function [Automatic restart], take appropriate measures to make sure that, in case of a possible unmonitored restart, no hazard for persons or risk of damage to property can occur.

In the as-delivered condition, the [Automatic restart] function is deactivated and has to be activated or deactivated manually.

The evaluation of potential risks by a possible unmonitored automatic restart as well as of appropriate security measures is the sole responsibility of the user!

Operation

WARNING

Whenever possible, use closed vessels for processing corrosive, toxic or biohazardous substances and seal them safely.

Increase the speed only gradually to the desired shaking frequency, especially with open vessels, and observe the fluid movements.

Please note the maximum lifting capacity of the platform of your device.

Lower the shaking frequency or reduce the total load on the platform when the device starts to vibrate during operation.

Always ensure that the vessels are evenly distributed on the platform.

Pay particular attention to an even distribution of the weight load when vessels of different sizes and/or differently filled vessels are placed on the platform at the same time!

Before switching on the device, make sure that all vessels are properly fixed on the platform.

Use suitable attachments for the used vessels. Further information on available accessories can be found on our website at www.heidolph.com!

Always keep existing ventilation slots clear to prevent the appliance from overheating. Close existing protective covers to prevent liquids and/or foreign objects from entering through appliance openings. Always use the necessary and appropriate personal protective equipment!

Secure the device in remote mode with a clearly visible warning sign and, if necessary, take further ambient-specific protective measures that protect against damage to property and injuries in the event of unexpected/unmonitored starting of the device.

Device service

DANGER

Switch the device's main switch off and disconnect it from the power supply before carrying out maintenance work, cleaning, or repairs.

When cleaning, avoid the penetration of liquids.

Before replacing the fuses, switch off the device and disconnect the power supply.

Always replace the two device fuses in pairs with original manufacturer fuses. Further information on available accessories can be found on our website at www.heidolph.com!

After fuse replacement, check the device for a safe condition according to IEC 61010-1.

CAUTION

Clean the device's surfaces with a soft, lint-free and only slightly moistened cloth.

Never use any aggressive or abrasive cleaning agents or aids.

Repairs

Repairs to the device may always be carried out by authorized experts! Unauthorized repairs during the warranty period will result in the loss of the warranty claim. The owner is generally liable for damage caused by unauthorized repairs. If repairs are necessary, contact an authorized dealer or our technical service.

Maintenance

There are no user-serviceable components in the unit housing. If necessary (in the event of abnormal operating behavior such as excessive noise or heat generation, for example) contact our technical service.

Check the caps and fasteners of all glass components for correct and tight fit at regular intervals and especially before each start-up of the device.

Check the level sensor in the heating bath regularly for proper operation, contamination, and limescale build-up to prevent malfunctions.

CAUTION

At loose/leaky connections, ambient air is sucked in during the vacuum build-up and the required vacuum pressure cannot be reached in a stable manner!

In the event of excess pressure, there is a risk that liquid will escape from the system! The device and surrounding areas can be severely contaminated by leaking media!

Tighten loose caps/fasteners by hand or correct the seating of the retaining clips.

Disposal

When disposing of the device, observe the provisions of the WEEE Directive 2012/19/EU and its transposition into national law in the country of use.

When disposing of portable batteries, observe the provisions of the European Battery Directive 2013/56/EU and their implementation in national law in the country of use.

Check the device and all components for residues of substances that are hazardous to health, the environment and biohazardous before disposal.

Properly remove and dispose of residues of substances that are hazardous to health, the environment and biohazardous!

Contact details

Heidolph Scientific Products GmbH

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© Heidolph Scientific Products GmbH